# Operating Manual

Original instructions - January 2025

LSPone syringe pump

Advanced Microfluidics SA

LSPone

Copyright © 2016

Published by:  
ADVANCED MICROFLUIDICS SA  
Chemin de la Dent d'Oche 1A  
1024 Ecublens  
Switzerland

[www.amf.ch](http://www.amf.ch)

*Version 3.2, en*

The information and data contained in this documentation is subject to change without notice. The reproduction, distribution and utilization of this document as well as the communication of its contents to others without explicit authorization is prohibited.

All rights reserved in the event of the grant of a patent, utility model or design.

The general terms and conditions of *Advanced Microfluidics SA* apply. Alternative agreements must be in written form.

*Advanced Microfluidics SA* shall not be liable for errors contained in this document or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

# Contents

<b>1</b>	<b>Introduction</b> .....	<b>7</b>
<b>1.1</b>	<b>Foreword</b>	<b>7</b>
<b>1.2</b>	<b>Application Purpose</b>	<b>7</b>
<b>1.3</b>	<b>Regulatory Considerations</b>	<b>7</b>
1.3.1	Declaration of Conformity CE .....	8
1.3.2	CB Test Certificate .....	9
1.3.3	English - Regulatory Considerations .....	10
1.3.4	Français - Considérations réglementaires .....	11
1.3.5	Deutsch - Regulatorischen Betrachtungen .....	12
1.3.6	Español - Consideraciones regulatorias .....	13
1.3.7	Italiano - Considerazioni normative .....	14
1.3.8	Nederlands - Regelgevende Overweging .....	15
1.3.9	Português – Considerações reguladoras .....	16
1.3.10	PYCKKNÑ .....	17
<b>1.4</b>	<b>Safety Considerations</b>	<b>18</b>
1.4.1	English - Safety considerations .....	19
1.4.2	Français - Considérations de sécurité .....	21
1.4.3	Deutsch - Sicherheit Betrachtungen .....	23
1.4.4	Español - Consideraciones de seguridad .....	25
1.4.5	Italiano - Considerazioni di sicurezza .....	27
1.4.6	Nederlands - Veiligheids Overwegingen .....	29
1.4.7	Português – Considerações de segurança .....	31
1.4.8	PYCKKNÑ .....	33
<b>1.5</b>	<b>Functional Description</b>	<b>34</b>
1.5.1	General Description .....	34
1.5.2	Syringe and syringe drive .....	34
1.5.3	Valve module .....	35

## I Hardware


<b>2</b>	<b>Hardware description</b> .....	<b>38</b>
<b>2.1</b>	<b>Glossary</b>	<b>38</b>
<b>2.2</b>	<b>Cabling</b>	<b>38</b>
2.2.1	Power supply .....	38
2.2.2	Female Mini-USB .....	39
2.2.3	Female DB9 connector .....	39
<b>2.3</b>	<b>Chemical compatibility</b>	<b>40</b>
<b>3</b>	<b>Valve module description</b> .....	<b>41</b>
<b>3.1</b>	<b>Low and small carryover volume valves</b>	<b>41</b>
3.1.1	Description .....	41
3.1.2	Port numbering .....	41
<b>3.2</b>	<b>Ultra low carryover volume valves</b>	<b>42</b>
3.2.1	Description .....	42
3.2.2	Port numbering .....	43
3.2.3	Standard use of the 6-port ultra-low carryover volume valve .....	44
<b>3.3</b>	<b>Minimal carryover volume operation</b>	<b>45</b>
3.3.1	Carryover volume .....	45
3.3.2	Simplified priming process .....	45
<b>4</b>	<b>Maintenance operations</b> .....	<b>47</b>
<b>4.1</b>	<b>Good practices</b>	<b>47</b>
<b>4.2</b>	<b>Cleaning procedure</b>	<b>47</b>
<b>4.3</b>	<b>Gas bubbles</b>	<b>49</b>
4.3.1	Bubble sources .....	49
4.3.2	How to avoid/reduce the creation of nuclei .....	50
4.3.3	Remove the trapped bubbles .....	51
<b>4.4</b>	<b>Changing the syringe</b>	<b>52</b>
<b>5</b>	<b>Hardware configurations</b> .....	<b>55</b>
<b>5.1</b>	<b>Valve configuration</b>	<b>55</b>
<b>5.2</b>	<b>Syringe configuration</b>	<b>55</b>

## II Software


<b>6</b>	<b>Software Communication</b> .....	<b>58</b>
<b>6.1</b>	<b>Communication Protocol</b>	<b>58</b>
6.1.1	Data Terminal Protocol .....	58
6.1.2	Using DT Protocol with Microsoft Windows .....	58
6.1.3	LSPone Command Set .....	60

---

<b>6.2</b>	<b>Debugging software problems</b>	<b>68</b>
6.2.1	Visual feedback to the user . . . . .	68
6.2.2	Error Codes and Pump Status . . . . .	69
6.2.3	Troubleshooting . . . . .	73
<b>7</b>	<b>GUI LSPoneQuick . . . . .</b>	<b>74</b>
7.1	Description	74
7.2	Requirements	74
<b>8</b>	<b>Labview VI . . . . .</b>	<b>75</b>
8.1	Description	75
8.2	Requirements	76
<b>9</b>	<b>Python scripting . . . . .</b>	<b>77</b>
9.1	Description	77
9.2	Requirements	77
<b>A</b>	<b>Resolution . . . . .</b>	<b>78</b>
A.1	Definitions	78
A.2	Flow rate in function of plunger speed	79
A.3	Volume in function of plunger displacement	79

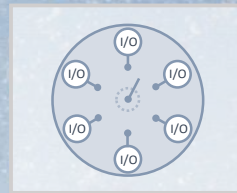# 1. Introduction
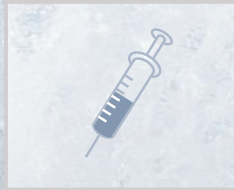
## 1.1 Foreword

The *Advanced Microfluidics SA*'s team thanks you for acquiring a LSPone syringe pump. This operating manual is dedicated to help you make the most of this device and thus bring success to your microfluidic projects. We remain available for any questions or suggestions that you may have.

The LSPone syringe pump system may only be operated after carefully reading and understanding this manual.

## 1.2 Application Purpose

The LSPone syringe pump is a programmable, precision liquid handling system. It is designed to enable pipetting, diluting, and dispensing of different liquids. It is controlled by an external computer or microprocessor and is powered by a 18-24V power supply. This pump is usually used in a laboratory-like environment

## 1.3 Regulatory Considerations

## EC Declaration of Conformity

According to EC directive 2006/42/EC on machinery (Annex II A)

The equipment which accompanies this declaration is in conformity with EU Directive(s):-  
*2006/42/EC Machinery Directive*  
*2006/95/EC Low Voltage Directive*  
*2004/108/EC Electromagnetic Compatibility Directive*

This declaration relates exclusively to the machinery in the state in which it was placed on the market, and excludes components which are added and/or operations carried out subsequently by the final user. The declaration is no more valid if the product is modified without agreement

### Manufacturer:

Advanced Microfluidics SA  
Chemin de la Dent d'Oche 1A  
1024 Ecublens  
Switzerland

A copy of the Technical file for this equipment is available from:

Remy Rysman  
Advanced Microfluidics SA  
Chemin de la Dent d'Oche 1A  
1024 Ecublens  
Switzerland

### Description of Equipment:

LSP one Syringe Pump

Model/type:

LSP ONE

The following harmonized standards have been used:-

*EN 61326-1:2013, IEC 61326-1:2012 (ed2.0)*

*EN 61010-1*

*EN ISO 14121-1*

Only for EN 61326-1:2013, IEC 61326-1:2012 (ed2.0):

*A sample of this machinery has been presented to Notified Body number STS 0024.*

*Electrosuisse, Montena EMC, route de Montena 75, 1728 Rossens, who have issued an EC type-examination certificate Number **16-MO-0065.E01** dated 2016.11.23.*

*The equipment in respect of which this declaration is made conforms to the example to which that certificate relates, and that certificate remains valid.*

Authorized signatory of manufacturer:

Ecublens, 23.11.2016

Barraud, Antoine / CEO

Place, Date

Surname, First name and function of the  
signatory

Signature

IEC SYSTEM FOR MUTUAL RECOGNITION OF TEST CERTIFICATES FOR ELECTRICAL EQUIPMENT (IECEE) CB SCHEME

**CB TEST CERTIFICATE**

Product	Syringe pump for academic research laboratories		
Name and address of the applicant	Advanced Microfluidics SA	Chemin de la Dent d'Oche 1a 1024 Ecublens VD	SWITZERLAND
Name and address of the manufacturer	Advanced Microfluidics SA	Chemin de la Dent d'Oche 1a 1024 Ecublens VD	SWITZERLAND
Name and address of the factory	Advanced Microfluidics SA	Chemin de la Dent d'Oche 1a 1024 Ecublens VD	SWITZERLAND
<i>Note: When more than one factory, please report on page 2</i>	<input type="checkbox"/> Additional Information on page 2		
Ratings and principal characteristics	LSP One: 18 VDC  AC/DC Switching Adaptor MEAN WELL (GSM40A18-P1J): Input: 100-240 VAC., 50/60 Hz, 1.0-0.05 A Output: 18 VDC, 2.22 A, 40 W max. Class I, IP20		
Trade mark (if any)	Advanced Microfluidics		
Customers's Testing Facility (CTF) Stage used	---		
Model / Type Ref.	LSP One GSM40A18-P1J		
Additional information (if necessary may also be reported on page 2)	---		
	<input type="checkbox"/> Additional Information on page 2		
A sample of product was tested and found to be in conformity with IEC	IEC 61326-1:2012		
National differences	EU Group Differences; EU Special National Conditions; EU A-Deviations; AU; CA; ID; IL; JP; KR; MY; SG; US; ZA		
As shown in the Test Report Ref. No. which forms part of this Certificate	16-MO-0065.E01 + .E02		

This CB Test Certificate is issued by the National Certification Body

Electrosuisse  
Luppmenstrasse 1  
8320 Fehraltorf  
SWITZERLAND

Signed by: Martin Plüss  
Date: 2016-11-30

### 1.3.3 English - Regulatory Considerations

#### Description

The LSPone is a general laboratory module. It is not intended for medical purposes therefore not subject to FDA regulatory approval.

*The LSPone must not be used as a medical device or for medical purposes.*

#### CE

CE certification is valid for the LSPone product. This does not apply to prototypes that are lent for evaluation.

#### Equipment ratings

<b>Environment</b>	
Use & Operation	Indoor
Operating temperature	15 to 40°C (41 to 104°F)
Storage temperature	-20 to 70°C (-4 to 158°F)
Humidity range	20 to 80% non condensing
Altitude	0 to 2000 m
<b>Mechanical Data</b>	
Weight	2.2 kg
Dimensions (L x W x H)	143 x 85 x 245 mm
Shipping weight	3.4 kg
Shipping dimensions (L x W x H)	413 x 301 x 125 mm
<b>Electrical Data</b>	
Power requirements	90 to 260 VAC, 1A max, 50/60 Hz
Power rating	18-24 VDC, 2.2A, 40 W
<b>Interfaces</b>	
USB (mini connector)	1.1, 2.0 and 3.0
Serial	RS232 and RS485

TABLE 1.1 – Equipment ratings

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## 1.4 Safety Considerations

### 1.4.1 English - Safety considerations

For proper handling and care of the LSPone, it is essential that the operating personnel follow the general safety procedures and safety instructions described in this manual.

#### Environment of use

- The LSPone syringe pump should only be used within the limits given by the equipment ratings
- The LSPone syringe pump is intended for indoor use only
- The LSPone syringe pump should not be used in an explosive atmosphere or with potentially explosive substances
- Transportation, storage or operation of the devices below 0°C with water in the fluid passages may cause damage to the modules
- Always work in a clean and open area to ease manipulation and avoid risk of injury

#### Intended use & operation

The LSPone syringe pump is designed for precise pipetting, diluting and dispensing operations in the 0.1  $\mu\text{L}$  to 1 mL range. Any other use is considered improper and may result in damage to the pump and/or unreliable test results.

Only the original approved parts and accessories may be used with the LSPone syringe pump. Any alterations or modifications to the instrument may be dangerous and will void the warranty.

#### Operating the LSPone

- When using the LSPone, Good Laboratory Practices (GLP) should be observed
- Users should wear protective clothing, safety glasses and protective gloves, especially if working with radioactive, biohazardous or harsh chemicals
- During the operation of the LSPone syringe pump, stand clear of moving parts
- Do not block moving parts
- Never try to remove valves, syringes or tubings when the pump or the valve is moving
- Never move the LSPone while it is in operation

#### Electrical safety

- The LSPone syringe pump must be disconnected from the power source when removing or replacing any mechanical components
- Use the provided power supply with the LSPone syringe pump (GSM40A18-P1J)
- Never use a damaged power cord
- Do not try to open the housing
- Do not power a device with a damaged housing

#### Precautions with radioactive, biohazardous or harsh chemicals

Please use great care when manipulating any chemical that is potentially damaging to your health.

*The LSPone syringe pump does not provide any user protection against radioactive, biohazardous or harsh chemicals*

- Always wear protective clothing, safety glasses and protective gloves
- Clean the fluid passages (including valve and syringe) immediately after using radioactive, biohazardous or harsh chemicals
- Make sure that the fittings are correctly tightened to avoid undesired leakage
- Make sure that the glass syringe is not damaged and correctly tightened

*Any liquid output can produce a squirt of liquid if the pump is programmed to push liquid through this very output at a high flow rate. Do not forget to protect yourself.*

Should an accidental spill occur, turn off the instrument and wipe it down with the appropriate disinfectant or chemical. Remember to take into account the nature of the spill and the necessary safety precautions.

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## 1.5 Functional Description

### 1.5.1 General Description

The device is made of the following parts:

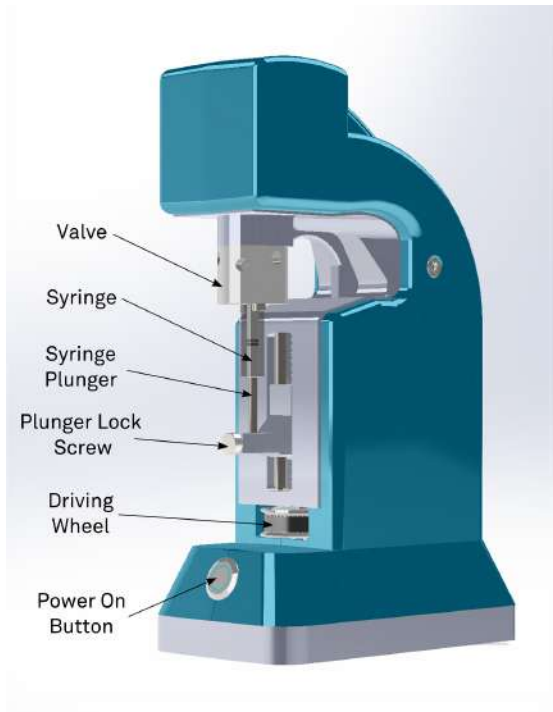
FIGURE 1.1 – LSPone syringe pump terminology 1/2

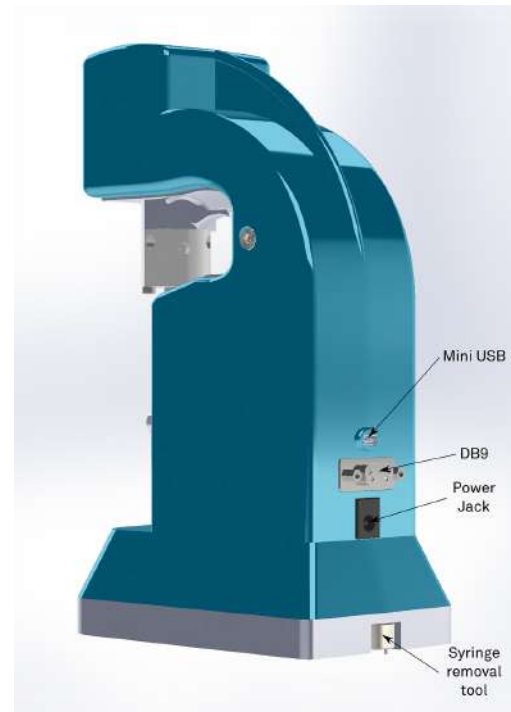
FIGURE 1.2 – LSPone syringe pump terminology 2/2

### 1.5.2 Syringe and syringe drive

The syringe assembly is made of a UHMW-PE or PTFE plunger tip which travels inside a glass syringe. The plunger is moved by a  $1.8^\circ$  stepper motor through an elastic transmission which enables a smoother displacement. The syringe plunger is tightened to the sliding part using a plunger lock screw.

The rated pressure for the syringe is 5 bars (0.5 MPa or 72.5 PSI).

FIGURE 1.3 – Detail of plunger fixation when correctly tightened

*Before performing any movements with the LSPone pump, please check that the plunger is correctly fixed to the slider. The screw should be completely inserted in the thread.*

The syringe volume is variable, between 25 and 1'000  $\mu\text{L}$ . For all syringes, the travel length of the plunger is 30 mm. When the pump is not powered, the syringe is removable. The plunger can also be displaced using the driving wheel only if the pump is not moving.

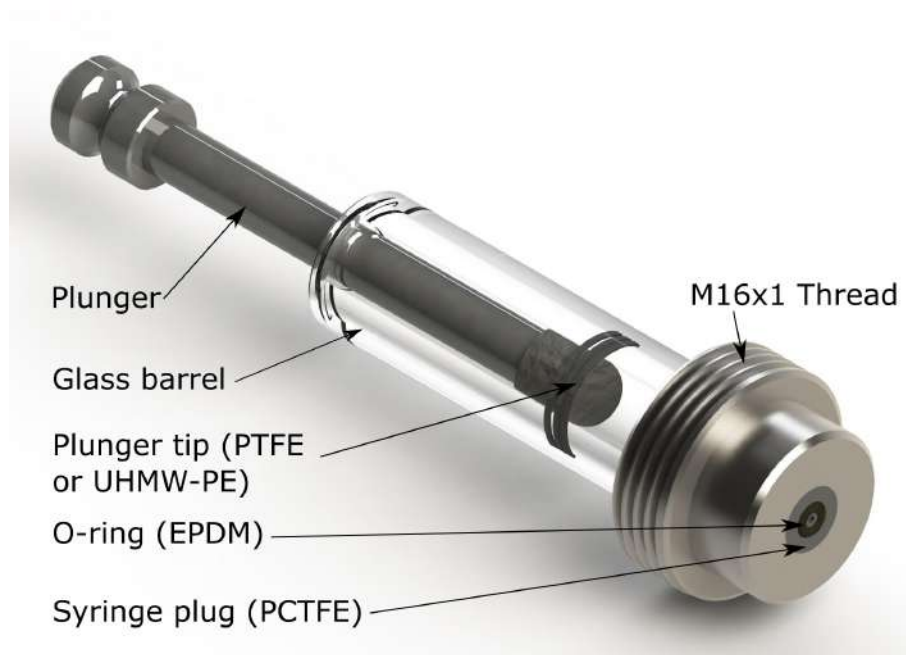
FIGURE 1.4 – Nomenclature for the syringe module

The PCTFE plug and the EPDM O-ring form the first and second sealing parts for the syringe. Therefore they should be treated with care. The O-ring should not be removed.

### 1.5.3 Valve module

The valve module enables the selection between the microfluidic ports (only one port is connected to the syringe at a given time). The valve module is made of fluoropolymers which have an outstanding chemical resistance. The plug which selects the connected channel is made of PTFE and the outer part, called valve seat, where the user can connect standard microfluidic fittings, is made of PCTFE.

The valve module, which is a closed loop system, works as follows:

- The user sends a command to connect a specific port
- The plug is rotated by a  $1.8^\circ$  stepper motor through the plug drive
- A position sensor detects very precisely the position of the plug thanks to indexes integrated in the plug itself
- Using this feedback the stepper motor can be stopped when the plug has reached the desired position

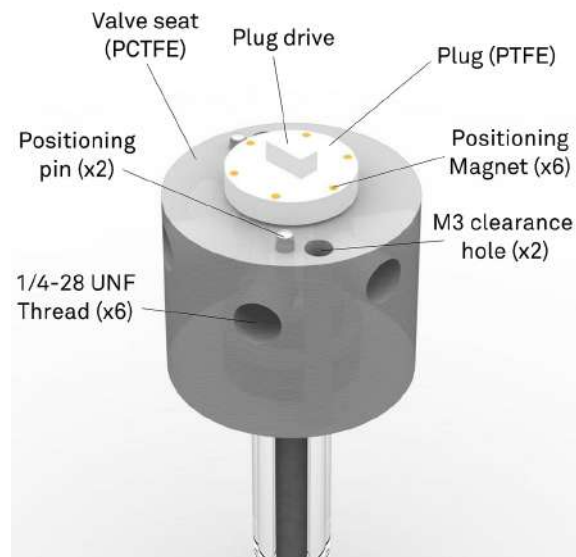
FIGURE 1.5 – Nomenclature for the valve module



When the pump is not powered, the valve module is removable by unscrewing the two M3x30 screws. The entire procedure is described in the maintenance section. The rated pressure for the valve module is 5 bars (0.5 MPa or 72.5 PSI).

# Hardware

<b>2</b>	<b>Hardware description</b> .....	<b>38</b>
2.1	Glossary	
2.2	Cabling	
2.3	Chemical compatibility	
<b>3</b>	<b>Valve module description</b> .....	<b>41</b>
3.1	Low and small carryover volume valves	
3.2	Ultra low carryover volume valves	
3.3	Minimal carryover volume operation	
<b>4</b>	<b>Maintenance operations</b> .....	<b>47</b>
4.1	Good practices	
4.2	Cleaning procedure	
4.3	Gas bubbles	
4.4	Changing the syringe	
<b>5</b>	<b>Hardware configurations</b> .....	<b>55</b>
5.1	Valve configuration	
5.2	Syringe configuration	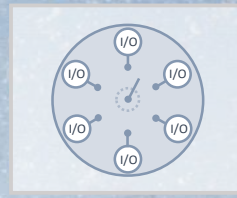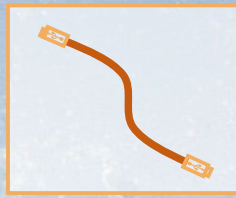

## 2. Hardware description

### 2.1 Glossary

**Carryover:** Contamination of a volume of fluid by residual fluid from a previous pick up or dispense. Carryover causes variability in the final volume and concentration.

**Carryover volume:** Volume of liquid that will be mixed with the next liquid. It is not stuck, but will be swept next time a liquid passes.

**Cavitation:** Formation of vapor cavities in a liquid – i.e. small liquid-free zones ("bubbles" or "voids") – that are the consequence of forces acting upon the liquid. It usually occurs when a liquid is subjected to rapid changes of pressure that cause the formation of cavities where the pressure is relatively low.

**Dead volume:** Volume that is “stuck” in the system (dead end), which is not cleanly swept and relies on diffusion to clear out.

**Internal volume:** Volume inside the system, from entrance to exit.

**Port:** The connection or orifice through which sample may be injected or dispensed.

**Priming:** Completely filling the pump tubing and syringe with bubble-free fluid to allow sustained, reproducible pumping action.

**Void volume:** Sum of the carryover volume and dead volume.

**Wetted surfaces:** The surfaces which are contacted by the sample stream.

### 2.2 Cabling

#### 2.2.1 Power supply

The LSPone syringe pump must be supplied with the AC/DC adapter which is provided with the pump. Using a different power supply is at the user's own risk.

FIGURE 2.1 – GSM40A18-P1J adapter from MeanWell provided with the pump

FIGURE 2.2 – Ferrite 600 04 27 from Sibalco to be added with 4 loops around the power cable

### 2.2.2 Female Mini-USB

The LSPone syringe pump can be controlled using a Mini-USB interface. In this case, the Mini-USB female connector works as a virtual serial port. The connection can be done by any mini-USB/USB cable provided that the provided ferrite of type 74271111S is locked around the cable.

FIGURE 2.3 – Ferrite 74271111S from Würth Elektronik to be locked around the USB cable connected to the LSPone

### 2.2.3 Female DB9 connector

The LSPone syringe pump can also be controlled using the DB9 connector<sup>1</sup>.

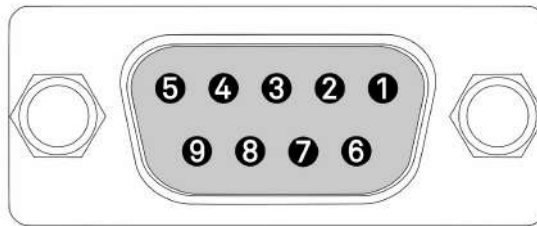
FIGURE 2.4 – DB9 pin-out on device side (LSPone)

Different communication protocols can be implemented with this single connector: RS-232, RS-485 or RS-422. Only the first two are implemented in the LSPone's software, but RS-422 can be implemented upon request.

Pin	Abbreviation	Name	Typical Purpose
1			
2	TX	Transmitted Data	Carries data from DCE (LSPone) to DTE
3	RX	Received Data	Carries data from DTE to DCE ( LSPone)
4	DTR	Data Set Ready	Not implemented. Connect to DSR
5	GND	Ground	Zero voltage reference
6	DSR	Data Terminal Ready	Not implemented. Connect to DTR
7	CTS	Clear To Send	Not implemented
8	RTS	Request To Send	Not implemented
9			

TABLE 2.1 – Pin assignment of the DB9 connector female on the LSPone for RS232 protocol

<sup>1</sup>Proper designation is DE-9 but the DB9 appellation is the most commonly used.

Pin	Abbreviation	Name	Typical Purpose
1			
2	D-	Data	Data. Negative driver
3			
4			
5	GND	Ground	Zero voltage reference
6			
7			
8	D+	Data	Data. Positive driver
9			

TABLE 2.2 – Pin assignment of the DB9 connector female on the LSPone for RS485 protocol

Pin	Abbreviation	Name	Typical Purpose
1			Not connected
2	TX-	Transmitted Data	Carries data from DCE (LSPone) to DTE. Negative output
3	RX+	Received Data	Carries data from DTE to DCE ( LSPone). Positive input
4			
5	GND	Ground	Zero voltage reference
6			
7	RX-	Received Data	Carries data from DTE to DCE ( LSPone). Negative input
8	TX+	Transmitted Data	Carries data from DCE (LSPone) to DTE. Positive output
9			Not connected

TABLE 2.3 – Pin assignment of the DB9 connector female on the LSPone for RS422 protocol

The pin assignment is always given from the point of view of the AMF module.

## 2.3 Chemical compatibility


The following parts can be in contact with the liquid:

1. The valve seat: PCTFE
2. The valve plug: PTFE or UHMW-PE
3. The syringe O-ring: EPDM
4. The syringe barrel: borosilicate glass
5. The syringe plunger tip: UHMW-PE or PTFE

These materials offer an excellent chemical compatibility. However Advanced Microfluidics SA cannot be held responsible for damages caused by the use of a specific chemical. Before using the device with a given chemical, please make sure that it is compatible with the aforementioned materials.

*The pump should not be used to process explosive or highly flammable fluids.*

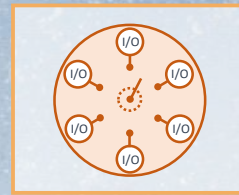## 3. Valve module description

The LSPone syringe pump can be used with several valve head types:

- Ultra-low carryover volume. Variants: 6 ports
  - Low carryover volume. Variants: 6, 8, 10, 12 ports
  - Small carryover volume. Variants: 8, 10 ports
- All of these valves have the particularity of having no dead volume.

### 3.1 Low and small carryover volume valves

#### 3.1.1 Description

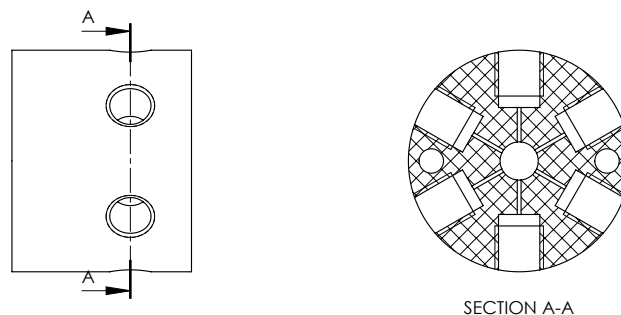
FIGURE 3.1 – Section view of the 6-port PCTFE valve seat revealing the microchannels

The 6-port low carryover volume valve (V-D-1-6-050-C-P) features 6 microfluidics ports that can be connected **one at a time** to the syringe content. The channels have a diameter of 0.5 mm in the valve seat and in the plug. Standard 1/4-28 UNF fittings can be connected to these channels.

The very small dimension of the channels inside the valve module reduces the carryover volume of the pump. The ports are organised on 1 stage, as shown in Figure 3.1, separated by 60°.

The same concept can be applied for the other models.

The internal and carryover volume increase for the 10 and 12-port valves, as they required a physically bigger valve head (otherwise we cannot fit that many ports around the circumference of the valve).

#### 3.1.2 Port numbering

You will find the port numbers for the VD1-6-050 in Figure 3.2 and Figure 3.3:

FIGURE 3.2 – Cut section revealing port numbering - VD1-6

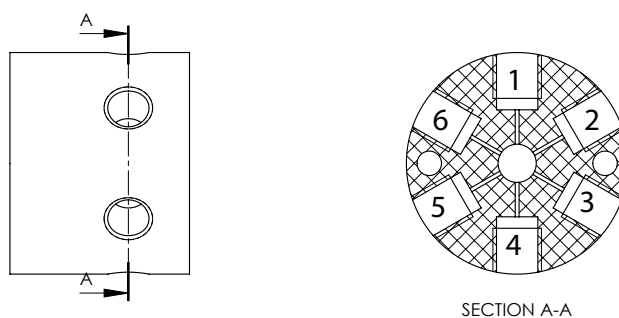
FIGURE 3.3 – Section view of the PCTFE valve seat with port numbering - VD1-6

**Note — Direction of rotation.**

The command I<n> corresponds to a clockwise rotation of 60° when incrementing the port number by 1. The command O<n> corresponds to a counterclockwise rotation of 60° when decrementing the port number by 1.

If you are on port 3, the command I<4> will be executed with a 60° rotation of the plug, while the command O<4> will be executed with a 300° rotation of the plug.

It is also possible to go to a port using the shortest way, regardless of the direction of rotation by using the command B<n>.

## 3.2 Ultra low carryover volume valves

### 3.2.1 Description

This very specific valve design has been patented for allowing minimal carryover volume. Depending on the requirements of the application, the valve can be used as a standard 6-port valve leading to a carryover volume of 3  $\mu\text{L}$  for the entire pump (in the worst case) or with a specific port choice for input and output, the pump can feature a carryover volume of 1.5  $\mu\text{L}$ .

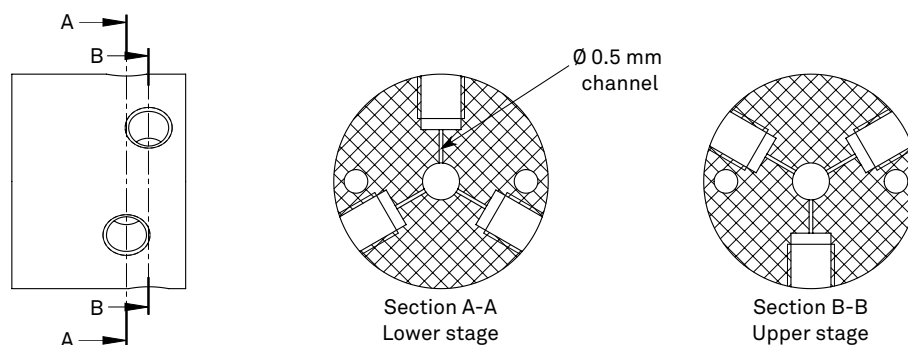
FIGURE 3.4 – Section view of the PCTFE valve seat revealing the microchannels - VD2-6

The 6-port ultra-low carryover volume valve features 6 microfluidics ports that can be connected **one at a time** to the syringe content. The channels have a diameter of 0.5 mm in the valve seat and in the plug. Standard 1/4-28 UNF fittings can be connected to these channels.

The very small dimension of the channels inside the valve module reduces the carryover volume of the pump. The ports are organized on 2 stages as shown in Figure 3.4. This configuration also minimizes the leakage rate since the ports are separated by 120°.

### 3.2.2 Port numbering

This feature requires a specific geometry for the microchannels and explains the specific port numbering depicted in Figure 3.5.

FIGURE 3.5 – Cut section revealing port numbering - VD2-6

Since the ports are located on 2 stages, the rotations from one port to the other when they are located on different stages may be counterintuitive. The port numbering is established as follows: between two consecutive ports (e.g.  $1 \Rightarrow 2$ ) the plug rotates in a clockwise manner of 60° (smallest possible rotation) unless asked to turn in the opposite direction (i.e. the counterclockwise direction). Additionally, odd-numbered ports are positioned below even-numbered ports according to Figure 3.6.

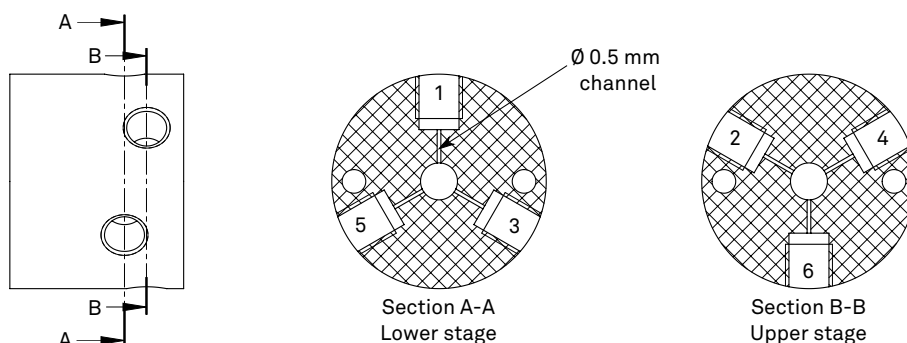
FIGURE 3.6 – Section view of the PCTFE valve seat with port numbering

**Note — Direction of rotation.**

The definition for rotating direction is given in Figure 3.7. The default rotation is triggered by the command `I<n>` which leads to a clockwise rotation of the plug. Moving from port 3 to port 4 will be done by a 60° clockwise rotation of the plug if called by the command `I<4>`. On the other hand, the user might want to rotate the plug in the counterclockwise direction with command `O<4>`. This will result in a 300° rotation of the plug for moving from port 3 to port 4.

FIGURE 3.7 – Direction of rotation definition for the plug - clockwise rotation corresponds to a 60° rotation when incrementing the port number by 1, counterclockwise rotation corresponds to a 60° rotation when decrementing the port number by 1

### 3.2.3 Standard use of the 6-port ultra-low carryover volume valve

In many applications, the valve can be used as a standard 6-port valve. Each port can be used as an input or an output similar to competitors' valves but the pump's carryover volume will remain below 5  $\mu\text{L}$ . This low carryover volume will prevent the appearance of an undesired air layer in the syringe.

In this mode, you can pick up up to 5 samples and dispense them through the remaining port for the desired microfluidics application.

**■ Application Notes**

You can find various application notes on our website at: <https://amf.ch/new/applications/>

### 3.3 Unleashing the minimal carryover volume capabilities of the ultra-low carryover volume valve

For some applications, the user will desire to use the specific feature of the LSPone syringe pump. The minimal carryover volume offered by our innovative construction enables maximum precision and reduced contamination between samples. Cleaning steps are thus made easier and more efficient.

#### 3.3.1 Carryover volume

The typical evidence of the carryover volume of a pump is the apparition of an air layer at the bottom of the plunger as illustrated in Figure 3.8. This air layer is tricky to remove and will cause inaccuracy in dispensed volumes. The LSPone syringe pump will already reduce this issue in the standard configuration since its carryover volume is extremely small. For more demanding applications, the pump offers a specific way to further reduce the carryover volume and its consequences.

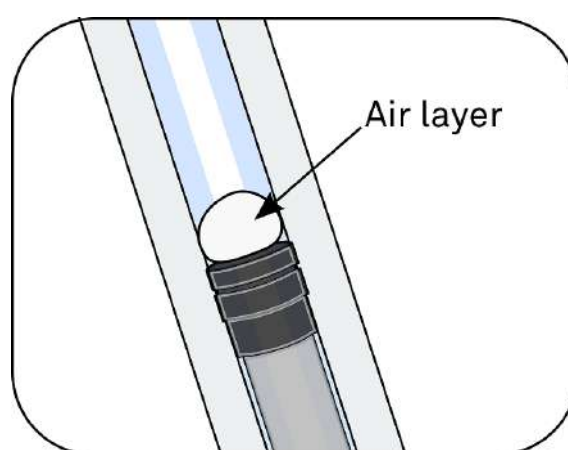
FIGURE 3.8 – Air layer appearing with conventional syringe pumps

#### 3.3.2 Simplified priming process

A minimal carryover volume offers several advantages. The chosen case, here, is the priming where the user wants to have the microfluidics tubings filled with the right reagent for his experiment. The priming step is usually an exhausting task with conventional syringe pumps due to the air layer.

The patented valve design offers a simplified procedure to prime your tubings before starting the experiment. This principle of operation is based on the two stages of the valve module. It should be performed with the input and the output ports on separate stages. The input tubing can be connected on the lower and the output port on the upper tubing or vice versa. The two solutions are equivalent. The complete process is described in Figure 3.9.


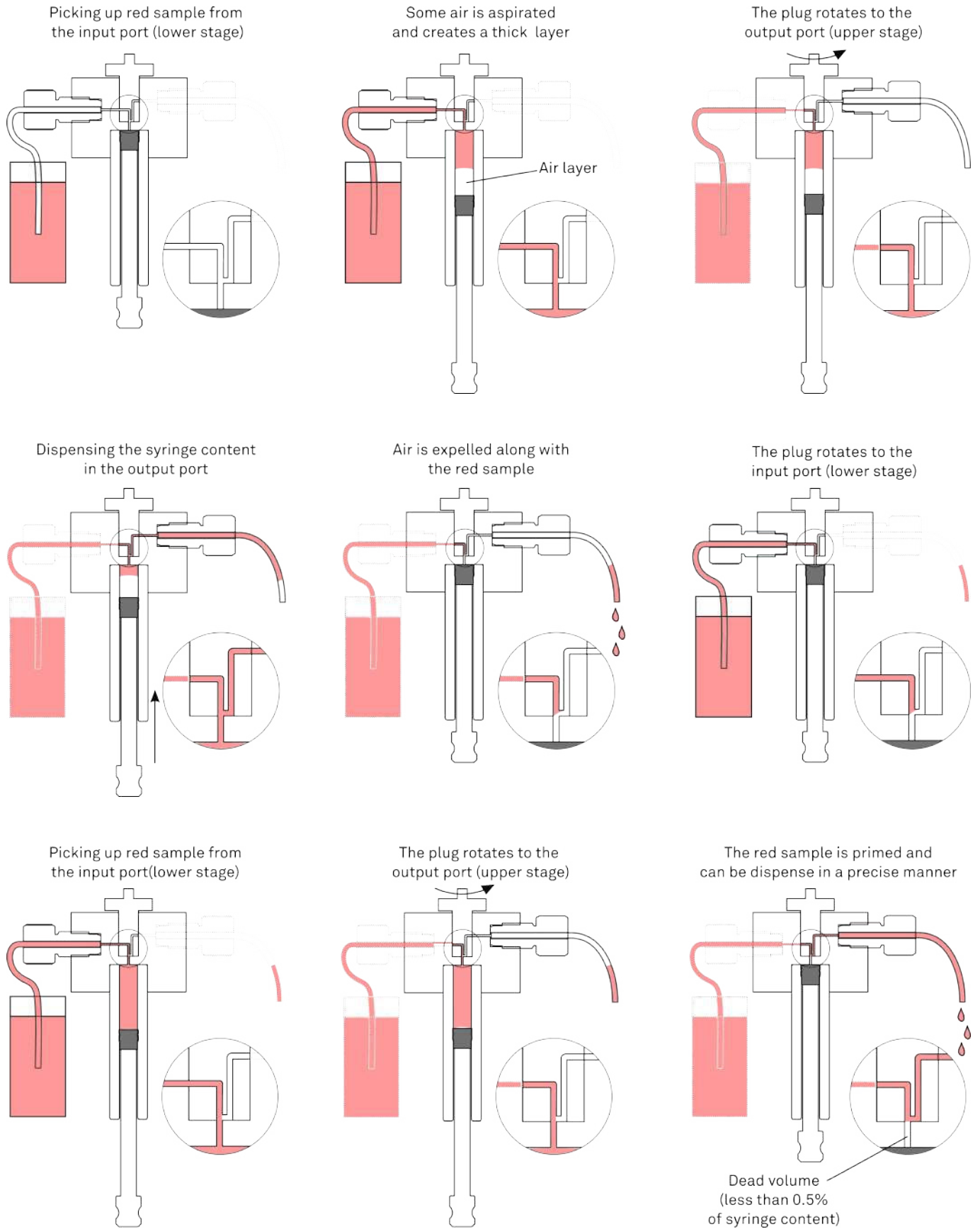
FIGURE 3.9 – Automated priming process with LSPone

## 4. Maintenance operations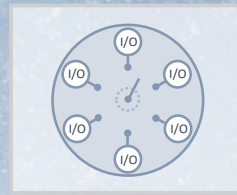

### 4.1 Good practices

In order to avoid premature damage of the LSPone or the valve, there is some good practice to have when you are using the pump:

- Try to avoid as much as possible dust entering the fluidic system because it can get block into the valve and create a groove in it. This would lead to leakage or bubble creation
  1. When your are screwing your tubing to the valve port, make sure to not have dust in the port. You can use dust free cleaning spray
  2. Filter your solutions before introducing them in the pump
  3. If you don't use one of the port of the valve, put a lid on it
  4. You can put the pump under a laminar flow to reduce the amount of dust
- Avoid over pressure, it can damage the valve or the syringe. Make sure the outlet is not blocked when you are dispensing
- Avoid solution that can crystallize in your microfluidic system. If you have to use this solution, clean your system after usage (see section 4.2)

### 4.2 Cleaning procedure

A cleaning procedure should be **performed at the end of each day or in between two different experiments**. An effective cleaning can be performed in the following manner:

- Bleach solution to kill microorganisms
  1. Prepare a solution of 1% chlorine bleach and deionized water
  2. Prepare a reservoir and place 1 tubing in the bleach solution and 5 tubings in a waste recipient
  3. Pick up the bleach solution through 1 port to entirely fill the syringe
  4. Sequentially output the syringe content through the remaining ports and throw it to the waste
  5. Repeat operation at least 2 times

**Note:** You can also use BD FACS clean solution that is typically used to clean flow cytometers
- Detergent solution to remove debris
  1. Prepare a solution of detergent (1% Mucosol) and deionized water
  2. Move the tubing from the bleach solution reservoir and place it into a reservoir of detergent solution
  3. Pick up the detergent solution through 1 port to entirely fill the syringe

4. Sequentially output the detergent solution through the remaining ports into the waste recipient
  5. Repeat operation at least 2 times
- Note:** 2% RBS 25 can also be used. For blood or plasma manipulation, you may use 1% Alconox. For the valve, you can make this washing step into a ultrasonic bath after a long period of using it.
- Rinsing water to prepare the pump for your next experiment
    1. Move the tubing from the detergent solution reservoir and place it into a reservoir of deionized water
    2. Pick up the rinsing solution through 1 port to entirely fill the syringe
    3. Sequentially output the rinse water through the remaining ports into the waste recipient
    4. Repeat operation at least 2 times
- Note:** Before the the rinsing step with the deionized water, you can add a washing step with a solvent such IPA or 70% ethanol (same manipulations as for the rinsing step)

**Note**

Thanks to our small internal volume, each cleaning cycle is extremely efficient.

## 4.3 Gas bubbles

Gas bubbles in a microfluidics setup is a very common issue and it can be quite difficult to address it. The bubbles can be trapped or directed into a waste output, but there is always a risk of some gas bubbles being created in the setup. The goal of the following section is to give you some hints to avoid them as much as possible and to remove them if they are trapped.

### 4.3.1 Bubble sources

#### Gas coming from leakage

The first possible source of gas bubble comes from the aspiration by the pumping system of external air. This is possible when some part of the microfluidics setup is not air tight. We propose a methodical approach to identify the leakage source. First check the following elements of the microfluidics setup:

- **Connectors:** Verify that every connector is tightly screwed/connected and that it suits well with the tubing you are using. We advise using connectors with a PTFE ferrule that is compressed when the connector is screwed. Make sure that the mating surfaces (PTFE ferrule and valve port surface) are clean and without scraps. To test if the leakage comes from a specific connector, spray a solution containing detergent around the connector and pressurize at 1 or 2 bars the connection with a syringe full of air for example. If there is a leakage, you will see some bubbles around the connector.
- **Valve:** Keep the solution you are passing through the valve as clean as possible. If dust or solid particles pass through your tubing and get stuck in the valve, it can create groove in the valve when this one is rotating. This will eventually lead to leakage and the valve will have to be changed. If you suspect that the valve is leaking, you can plug the port you are not using with a cap to see if it reduces the leakage. If the leakage is reduced, it means that the valve is damaged. After running a lot of cycles, the valve can wear out and the sealing between the rotor and the stator may be defective.
- **Syringe:** As for the valve, the syringe can wear out after a lot of cycles. Check the glass barrel for cracks especially after a choc. If you see a crack on the glass barrel, the syringe has to be replaced as the crack is a source of leakage. If the syringe looks good but you experience leakage, make sure that the small O-ring is correctly installed between the valve and the syringe.

As a general rule, to check if the valve or the syringe is the source of your bubble problem, change them one at a time and check if the situation is better.

#### Gas coming from the solution

The concentration of air molecule is given by the Henry's law:  $c_s = H(T)p_g$  where  $H(T)$  is the solubility constant (increase if the temperature  $T$  decrease) and  $p_g$  is the pressure of the gas in vapor phase. We have an equilibrium when the actual concentration of gas in the liquid  $c$  is equal to  $c_s$ . When the temperature or the pressure changes, there is some variation of  $c_s$  which leads to a non-equilibrium state described by:  $\zeta = c/c_s - 1$ . When  $\zeta < 0$ , we have a undersaturation situation and the gas wants to go in the solution (would reduce the size of the bubbles) and when  $\zeta > 0$ , we have a supersaturation situation and the gas wants to go out of the liquid and this makes bubbles growing. This supersaturated situation occurs when we reduce the pressure (by aspirating the liquid with a pump for example) or when we increase the temperature.

The bubbles created in the solution have a starting point called a nuclei and when this nuclei is formed, the bubbles start growing due to a supersaturated situation. Those nuclei can appear in the bulk of the solution if we have a strong supersaturated solution ( $\zeta \gg 0$ ) and this results in microbubbles flowing continuously through your tubing, or they can appear on the surface of your

tubing/connector/valve/syringe due to strong supersaturated solution (nuclei form on the surface and then in the bulk because the surface acts as a nuclei catalyzer) or due crevices or sharp edges during the wetting as shown in figure 4.1 (source : Pereiro, Iago, et al. "Nip the bubble in the bud: a guide to avoid gas nucleation in microfluidics." Lab on a Chip 19.14 (2019): 2296-2314).

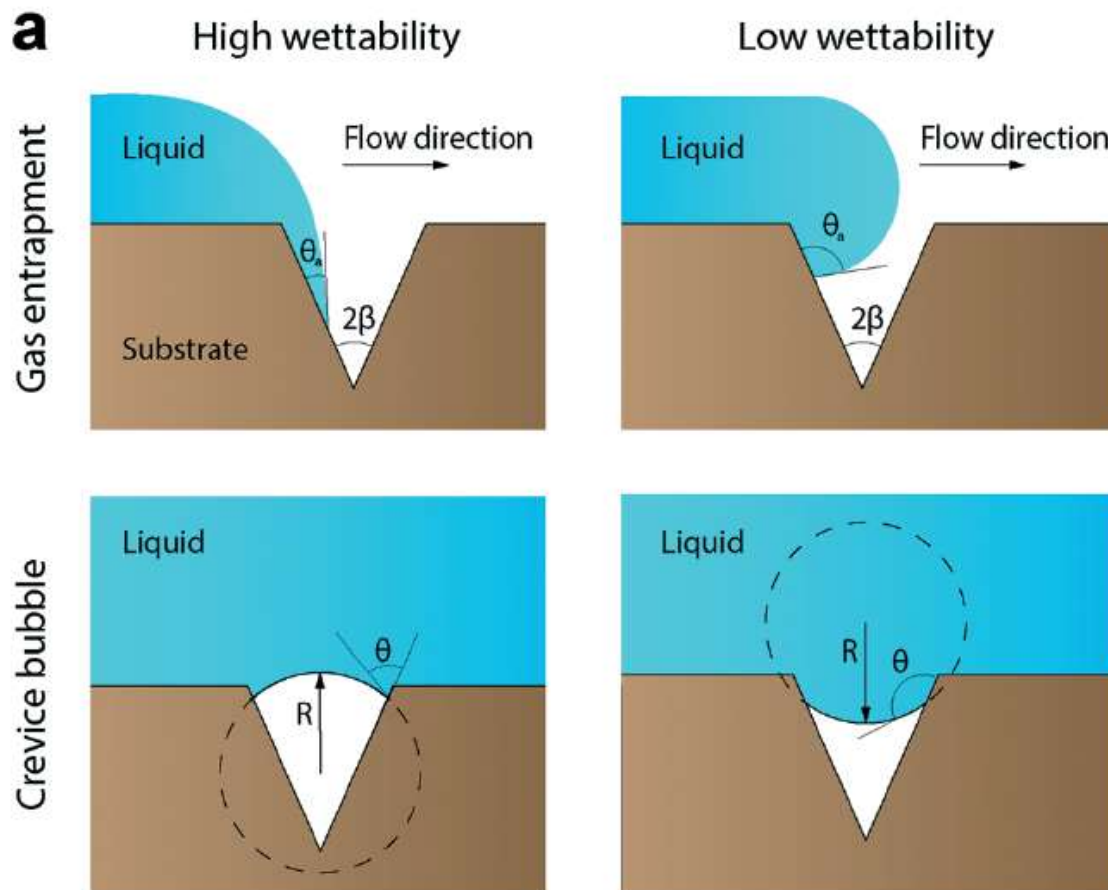
FIGURE 4.1 – Image that describe the formation of bubble due to a crevice during the wetting of the microfluidic network (source : Pereiro, Iago, et al. "Nip the bubble in the bud: a guide to avoid gas nucleation in microfluidics." Lab on a Chip 19.14 (2019): 2296-2314)

The crevice or sharp edge bubbles are the main source of "in solution created" bubbles. It acts as gas tank and when we have a supersaturation situation, the gas volume increases until the bubble is too big to stay trapped by the surface and part of bubble is released. This gas tank is a continuous source of bubble and more you are in a supersaturation situation, more often you will have bubble released from this gas tank. The surface nuclei are maintained on the surface due to the surface tension (the more hydrophobic the surface is, the more the gas bubble will stay trapped) and act as a trap for the microbubbles that are flowing through the tubing. Because of this we have an accumulation effect, the surface bubbles become bigger and bigger until they are too big to be maintained by the surface and they are then released.

#### 4.3.2 How to avoid/reduce the creation of nuclei

- Degas your solutions before using them and in the best case, keep them in this low gas concentration until they are used (for example, keep them under vacuum)
- Avoid big change of pressure and high velocity flow

- Try to keep your lab and your setup under a constant temperature (for example during the night, the temperature may decrease and re-increase in the morning leading to bubble creation)
- **Connectors:** Try to avoid as much as possible sharp edges and dead volume in your connectors and try to avoid change of diameter in your tubing
- **Valve:** As for the leakage, try to avoid dust and particles in the valve because it would create groove that are source of crevice nuclei and leakage. Also, try to fit the diameter of your valve port with the diameter of your tubing
- **syringe:** During the priming of your system, a bubble is created in the syringe due to the inner volume of the pump. This bubble may be washed away during the priming but can also remain trapped by the plunger (more hydrophobic than the syringe wall in glass), is never released and act as a bubble source under supersaturation situation. The solution for this is to prime your system with a solution that has a lower surface tension such as ethanol or SDS.
- **Wetting:** To avoid the crevice nuclei during the priming, it is important to prime your system with a low flow velocity. It also helps if you are using a solution with a lower surface tension such as ethanol or SDS for the wetting of your system

#### 4.3.3 Remove the trapped bubbles

To remove the bubbles trapped in your system, you can tap on the tubing or induce vibration to your system. The most efficient way we have found is to flow a large gas (air) bubble followed by a solution containing ethanol: The air train picks up all the trapped bubbles and the ethanol brings a good wetting on the surfaces.

## 4.4 Changing the syringe

The user has the possibility of changing the syringe of the LSPone syringe pump. Different volumes are offered and can be ordered to match the application's requirements. To extend the lifetime of the pump, please follow the instructions below carefully when changing the syringe.

The pump layout is given hereafter as a reminder, so as to simplify the syringe change.

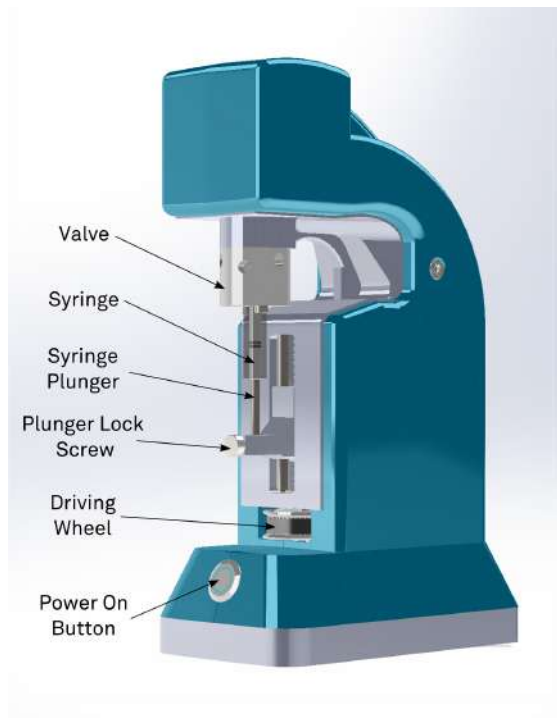
FIGURE 4.2 – LSPone syringe pump terminology 1/2

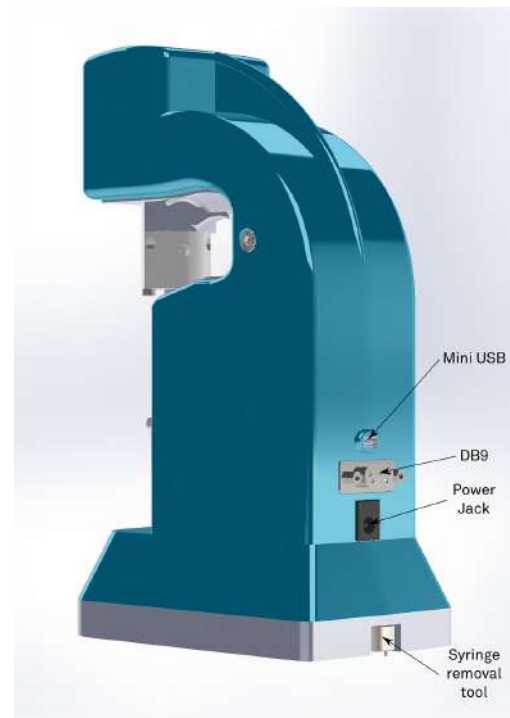
FIGURE 4.3 – LSPone syringe pump terminology 2/2

1. Disconnect the syringe pump from the power supply.
2. Position the syringe pump on an open and clean area.
3. Unscrew the plunger lock screw.
4. Use the driving wheel to be able to insert the syringe removal tool around the syringe barrel.
5. The pins of the syringe removal tool should be inserted in the two appropriate bores at the syringe basis. To remove the syringe unscrew the glass barrel using the removal tool. After many cycles, using pliers may be necessary (see Figure 4.4 and Figure 4.5)
6. After removal of the syringe, make sure that the black O-ring is correctly fitting in its groove and that it is not damaged.
7. Control the glass barrel. No crack or fissure should be seen on the glass.
8. If the removed syringe is in good condition, store it carefully as you will be able to use it again.
9. Perform the same quality check on the syringe that you plan to install.
10. Screw the syringe in the valve seat. Do not force the thread, the screwing process should be smooth.
11. Use the syringe removal tool to hand tighten the syringe. Do not use pliers or other tool when screwing the syringe.

FIGURE 4.4 – Insertion of the syringe removal tool around the glass barrel. Unscrew the syringe by turning in the highlighted direction.

FIGURE 4.5 – Syringe seal to be handled with care. Before screwing the syringe, ensure that the black O-ring is correctly inserted in the groove as shown here

*The syringe is designed to be tightened by hand. A good seal is only guaranteed if you follow this procedure.*

12. When the syringe is correctly tightened in position, use the driving wheel to come over the plunger.
13. Align the plunger bottom with the holder, according to Figure 4.6. The two surfaces marked in orange should be coplanar.
14. Screw the plunger lock screw and tighten it by hand.
15. Make sure that the plunger lock screw's thread is not visible anymore thus guaranteeing a safe fixation (see Figure 4.7 and Figure 4.8)

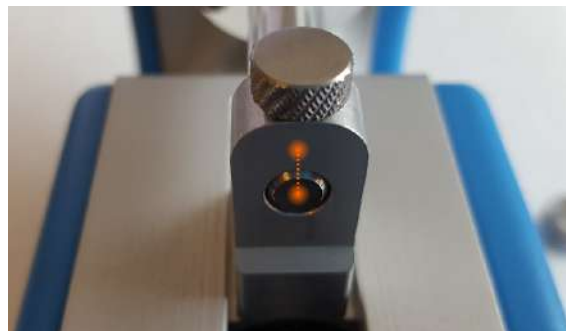
FIGURE 4.6 – Alignment to be obtained when screwing the plunger lock screw

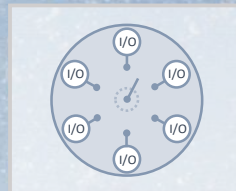

FIGURE 4.7 – Correct plunger lock screw fixation

FIGURE 4.8 – Incorrect plunger lock screw fixation

## 5. Hardware configurations
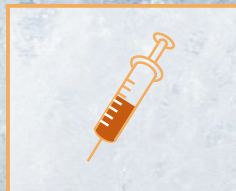
### 5.1 Valve configuration

Rated pressure: 5 bars

Reference	Nb of ports	Wetted materials	Internal vol.	Carryover vol.	Channel $\varnothing$
V-D-2-6-050-C-P	6	PCTFE, PTFE	5.2 $\mu\text{L}$	1.5 $\mu\text{L}$	0.5 mm
V-D-1-6-050-C-P	6	PCTFE, PTFE	3.5 $\mu\text{L}$	2.6 $\mu\text{L}$	0.5 mm
V-D-1-8-050-C-P	8	PCTFE, PTFE	3.5 $\mu\text{L}$	2.6 $\mu\text{L}$	0.5 mm
V-D-1-8-100-C-P	8	PCTFE, UHMW-PE	18.1 $\mu\text{L}$	11 $\mu\text{L}$	1 mm
V-D-1-10-050-C-P	10	PCTFE, UHMW-PE	4.5 $\mu\text{L}$	2.8 $\mu\text{L}$	0.5 mm
V-D-1-10-100-C-P	10	PCTFE, UHMW-PE	18.1 $\mu\text{L}$	11 $\mu\text{L}$	1 mm
V-D-1-12-050-C-P	12	PCTFE, UHMW-PE	4.5 $\mu\text{L}$	2.8 $\mu\text{L}$	0.5 mm

TABLE 5.1 – Valve types for LSPone

### 5.2 Syringe configuration

Rated pressure: 5 bars


Syringe Vol. [ $\mu\text{L}$ ]	Part Number	Plunger Material	Min. Flow Rate [ $\mu\text{L}/\text{min}$ ]	Max. Flow Rate [ $\mu\text{L}/\text{min}$ ]	Min Dosing volume [ $\mu\text{L}$ ]
25	S-25-P	PTFE	0.003725	750	0.05
50	S-50-P	PTFE	0.00745	1 500	0.1
100	S-100-P	PTFE	0.0149	3 000	0.2
100	S-100-U	UHMW-PE	0.0149	3 000	0.2
250	S-250-P	PTFE	0.03725	8 000	0.5
500	S-500-P	PTFE	0.0745	14 000	1
500	S-500-U	UHMW-PE	0.0745	14 000	1
1000	S-1000-P	PTFE	0.149	30 000	2

TABLE 5.2 – Syringe types for LSPone


# Software

<b>6</b>	<b>Software Communication</b> .....	<b>58</b>
6.1	Communication Protocol	
6.2	Debugging software problems	
<b>7</b>	<b>GUI LSPoneQuick</b> .....	<b>74</b>
7.1	Description	
7.2	Requirements	
<b>8</b>	<b>Labview VI</b> .....	<b>75</b>
8.1	Description	
8.2	Requirements	
<b>9</b>	<b>Python scripting</b> .....	<b>77</b>
9.1	Description	
9.2	Requirements	
<b>A</b>	<b>Resolution</b> .....	<b>78</b>
A.1	Definitions	
A.2	Flow rate in function of plunger speed	
A.3	Volume in function of plunger displacement	

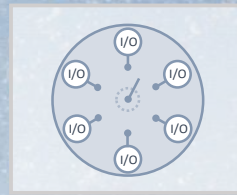## 6. Software Communication
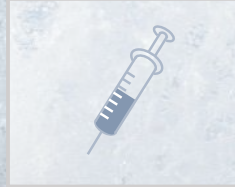
### 6.1 Communication Protocol

The LSPone syringe pump offers a data terminal protocol.

#### 6.1.1 Data Terminal Protocol

The data terminal protocol enables communication through the following pattern:

Character Format	
Parameter	Setting
Baud rate	9600
Data bits	8
Parity	None
Stop bit	1
Command Block	
1	Start command (ASCII “f”)
2	Pump address (1 by default, “_” is broadcast address)
3	Data block (length n)
...	
3+n-1	
3+n	Carriage Return (<CR>)
Answer Block	
1	Start answer (ASCII “f”)
2	Master address (ASCII “0”)
3	Error/Status character
4	Data block (length n)
...	
4+n-1	
4+n	ETX (0x03)
5+n	Carriage Return (<CR>)
6+n	Line feed (0x0A)

TABLE 6.1 – Data Terminal Protocol Format

Please note that due to the buffer size limitation, the length of the command block is limited to 512 characters. If necessary, the command string has to be split to avoid overflow. The looping functions will also help to reduce the number of required characters for a given sequence.

**Example 6.1** “/1ZR<CR>” will be sent to initialize the pump (see serial command description) In return, the pump will answer with “/0@<ETX><CR><LF>”. ■

#### 6.1.2 Using DT Protocol with Microsoft Windows

The LSPone can be controlled in the following manner:

1. Power on the pump (applying power will automatically switch on the pump).

2. Connect the LSPone to a communication port of the PC.
3. Open the *Device Manager* to identify which COM port is connected.
4. Open a Terminal or Terminal emulator such as RealTerm or HTerm.
5. Establish the connection by selecting the baud rate (9600), the correct communication port (here for example COM5), 8 data bits, 1 stop bit, no parity and no flow control.

FIGURE 6.1 – Settings for serial communication

6. Press *Change* to start the connection.
7. Send `/1ZR<CR>` with the *Send ASCII* button. The *End of Line* chars can usually be added automatically (here by ticking the `+CR` box).

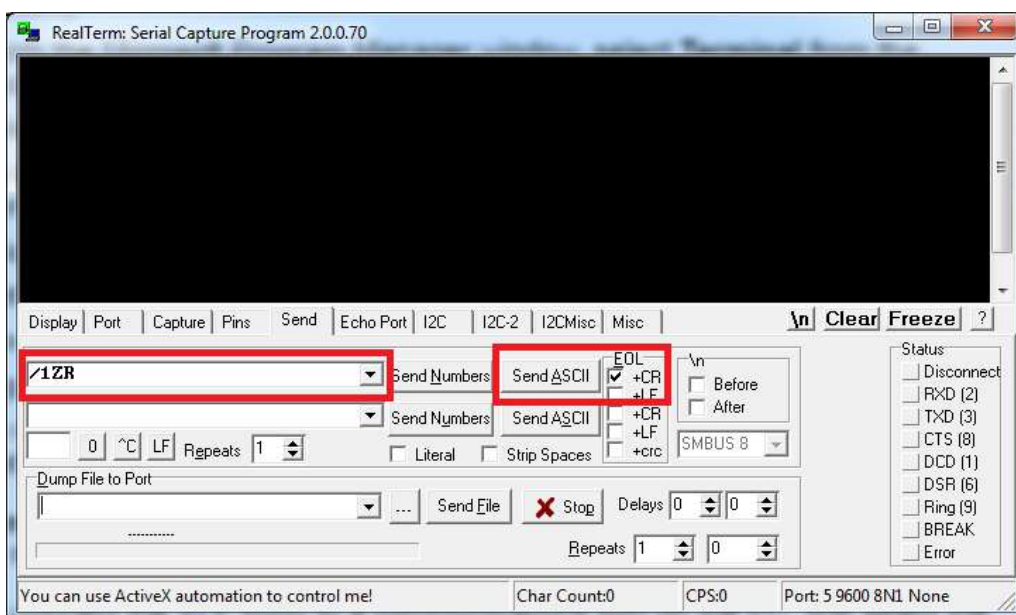
FIGURE 6.2 – Sending command configuration

8. Refer to the complete set of commands to use the pump according to your need.

### 6.1.3 LSPone Command Set

The LSPone simplifies your microfluidic automation thanks to a wide and well-tried command set. Many of the commands have default values; however, the default values may not provide the optimal settings for your application. Therefore it would be advantageous to spend some time exploring the many possibilities offered by the pump.

#### Command Execution Guidelines

- All commands, apart from those starting with "?" or "!" and from the abort command T, must be followed by the character [R] to be run.
- During the execution of a command, no new command is accepted, except for interrupt and report commands.
- The pump answers immediately after a command was sent. If an invalid command has been sent in the command string, the pump reports an error immediately.
- To reduce wear, avoid running the plunger dry in the glass cylinder.
- Keep your fingers out of the syringe slot for safety reasons.

#### Pump & Valve Configuration Commands

These commands are used to change your pump's setup. If you are unsure, please do not change these commands and ask us instead. You can also verify your setup using the report commands in 6.1.3.

Command	Operand Range	Default Operand	Power Up De-fault	Operand Description	Command Description
@ADDR=<n>	1..9 or A..E	1		1-character address	Pump address, "_" is broadcast address
!30<n>	0,1, 2, or 3	3		Plunger force	Setting of the plunger force "0" being high force, "1" being normal force, "2" being medium force, "3" being low force (no trailing [R])
!50<n>	0,1 or 2	2		Answer mode	Synchronous mode "0" : answer once when the command has been processed / Asynchronous mode "1" : answer when the command starts, for each query and when command is completed / Asynchronous mode + command counter "2" : same as asynchronous but add number of subcommand processed in its last answer (no trailing [R])
!80<n>	6, 8, 10 or 12	6		Number of valve positions, No trailing [R]	Valve configuration (no trailing [R])
!17					Reset valve motion counter (no trailing [R])
@RS232					Activate RS232 communication or serial-over-USB communication (activated by default) and deactivate RS485 communication
@RS485F					Activate RS485 communication (RS485 communication will not work if the mini USB cable remains plugged)

*We strongly advise using low force only when a small volume syringe (25, 50 or 100 µL) is used. Medium or normal forces can be used for syringe of 250 and 500*

*μL. High force is adapted to 1 mL-syringe and for applications exhibiting high fluidic resistance.*

When RS485 communication mode is activated, the module will not answer nor respond to query addressed via broadcast. But any command sent via broadcast will be performed. Hence if several modules are connected on a RS485 communication line, all the modules can be controlled with the broadcast address but no answer will be sent. In order to receive an answer from a specific module, each module has to have a unique address and it has to be addressed specifically.

When the module has RS485 communication activated, the firmware cannot be updated using the LSPoneQuick software.

*If a pump is not answering to requests sent by the broadcast command, it is likely that the module has RS485 communication activated.*

#### Example 6.2 How to use answer mode:

Answer mode	Description	Example	Answer	Timing
Synchronous	Only answer once when the command has been processed, can't have embedded queries	/_P100D50R	/0@	Immediate
Asynchronous	Answer when command starts, for each query and when command has ended, can have embedded queries	/_P100?4?49D50R	/0@ /0'100 /0c /0'	Immediate + after each query + after last command
		/_P100D50R	/0@ /0'	
Asynchronous + command counter	Same as asynchronous but add number of subcommand processed in its last answer	/_P100?4?49D50R	/0@ /0'100 /0c /0'4	Immediate + after each query + after last command
		/_P100D50R	/0@ /0'2	

## Control Commands

Command	Operand Range	Default Operand	Power Up Default	Operand Description	Command Description
R	N/A	N/A			Execute command string
X	N/A	N/A		No trailing [R]	Re-execute last executed command string
G<n>	0..60000	0		0 = Loop forever	Repeat command sequence
g	N/A	N/A		Loop depth = max 10	Mark the start of a repeat sequence
M<n>	0..86400000			Milliseconds	Delay command execution
H	N/A	N/A		No trailing [R]	Halt command - Pause the sequence after finishing the current move. The paused sequence will resume with "/IR<CR>" or alternatively the user can send a new command. The command H can also be integrated in a sequence to introduce a pause before the execution of the remaining commands
T	N/A	N/A		No trailing [R]	Hardstop - Interrupt the current move and suppress it from the sequence. The interrupted sequence will resume with "/IR<CR>" after the interrupted command or alternatively the user can send a new command
@POWEROFF	N/A	N/A			Shut down the pump

**Example 6.3** "/1gP2000D2000G3R<CR>" will trigger 3 back-and-forth 2000-step moves from the plunger.

"/1P2000HD2000R<CR>" will trigger a pick-up move of 2000 steps. The pump will then wait for "/IR<CR>" before executing the 2000-step dispense. The Halt command can also be executed on the fly. ■

## Initialization Commands

Command	Operand Range	Default Operand	Power Up Default	Operand Description	Command Description
Z<n>	(void) or 0..3	0		0 or (void)= high force, 1= normal force, 2=medium force, 3=low force	Initialize the plunger drive (with high, normal, medium or low plunger force) and home the valve
Y<n>	(void) or 0..3	0		0 or (void)= high force, 1= normal force, 2=medium force, 3=low force	Initialize the plunger drive (with high, normal, medium or low plunger force) and home the valve

*In our case, there is no difference between the two initialisation commands. Both exist to be compatible with the Cavro standard.*

## Valve Commands

Command	Operand Range	Default Operand	Power Up Default	Operand Description	Command Description
O<n>	1..6	N/A		Counter-clockwise plug movement (valve seen from above)	Move to valve port
I<n>	1..6	N/A		Clockwise plug movement (valve seen from above)	Move to valve port. Port numbering has been configured so that incrementing the argument of I by 1 triggers a plug rotation of 60 degrees only
B<n>	1..6	N/A		Shortest plug movement	Move to valve port with shortest path clockwise direction preferred

**Example 6.4** “/1M10000I2R<CR>” will be sent to wait 10 seconds before moving the plug to port 2 in a clockwise direction. In return, the pump will answer with “/0@<ETX><CR><LF>”

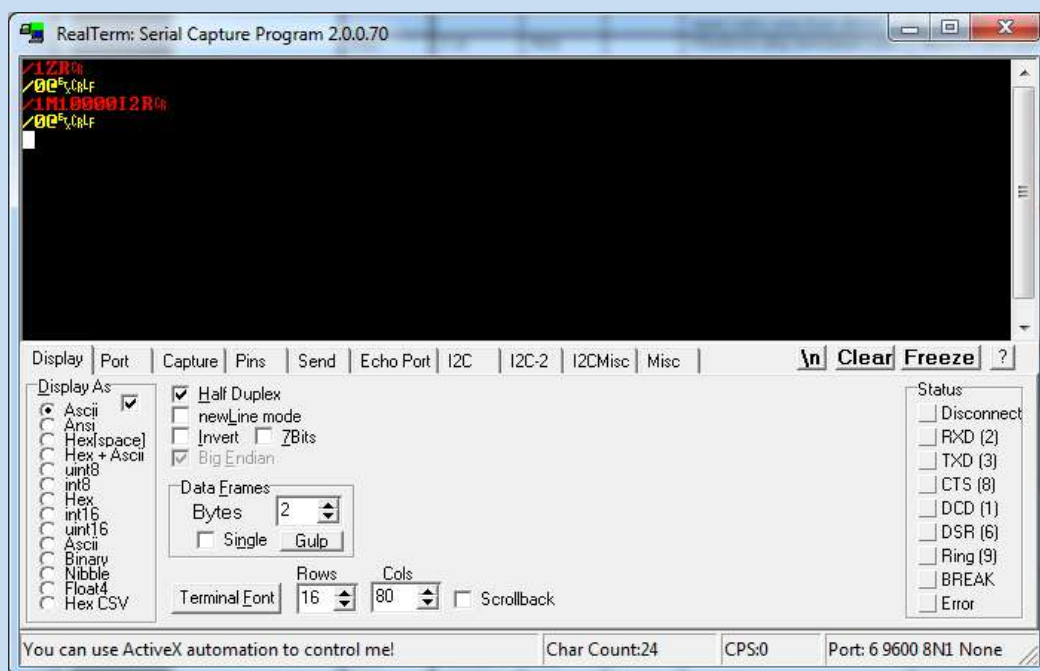
FIGURE 6.3 – Sending command via the terminal after starting the pump

Note in this example that the pump requires a homing command after power up.

## Plunger Movement Commands

Command	Operand Range	Default Operand	Power Up Default	Operand Description	Command Description
A<n> a<n>	0..3000 with N=0 0..24000 with N=1	N/A			Absolute Position
P<n> p<n>	0..3000 with N=0 0..24000 with N=1	N/A			Relative Pickup
D <n> d<n>	0..3000 with N=0 0..24000 with N=1	N/A			Relative Dispense

*In our case, there is no difference between the upper and lower caps for these positioning commands. Both exist to be compatible with the Cavro standard.*

## Set Commands

Command	Operand Range	Default Operand	Power Up Default	Operand Description	Command Description
L<n>	100..59590	N/A	1557	Acceleration rate (Pulses/sec <sup>2</sup> )	Set acceleration rate
l<n>	100..59590	N/A	59590	Deceleration rate (Pulses/sec <sup>2</sup> )	Set deceleration rate
N<n>	0..1	0		<0> = Microstep 0.01 mm resolution <1>= Microstep 0.00125 mm resolution	Scaling of dispense/pickup arguments. The motor is always driven in microstep mode
S<n>	10..40	N/A	22	See Table 6.2	Set peak speed code Only present to be compatible with other syringe pumps. Does not offer as many possibilities as the command V.
u<n>	1..214750	N/A		Peak speed (0.00745 Pulses/sec = 74.5 nm/s)	Set peak speed with ultra low flow rate resolution
U<n>	1..32000	N/A		Peak speed (0.05 Pulses/sec = 500 nm/s)	Set peak speed with low flow rate resolution Setting the speed with a factor 20 compared to V command
V<n>	0..1600	N/A	150	Peak speed (Pulses/sec = 0.01 mm/s)	Set peak speed V0 is a special setting and corresponds to 0.5 Pulse/sec

*The peak speed may not be reached depending on acceleration/deceleration settings and length of the movement*

**Example 6.5** “/1N1R<CR>” will be sent to choose the 0.00125 mm resolution microstepping mode. In return, the pump will answer with “/0‘<ETX><CR><LF>”

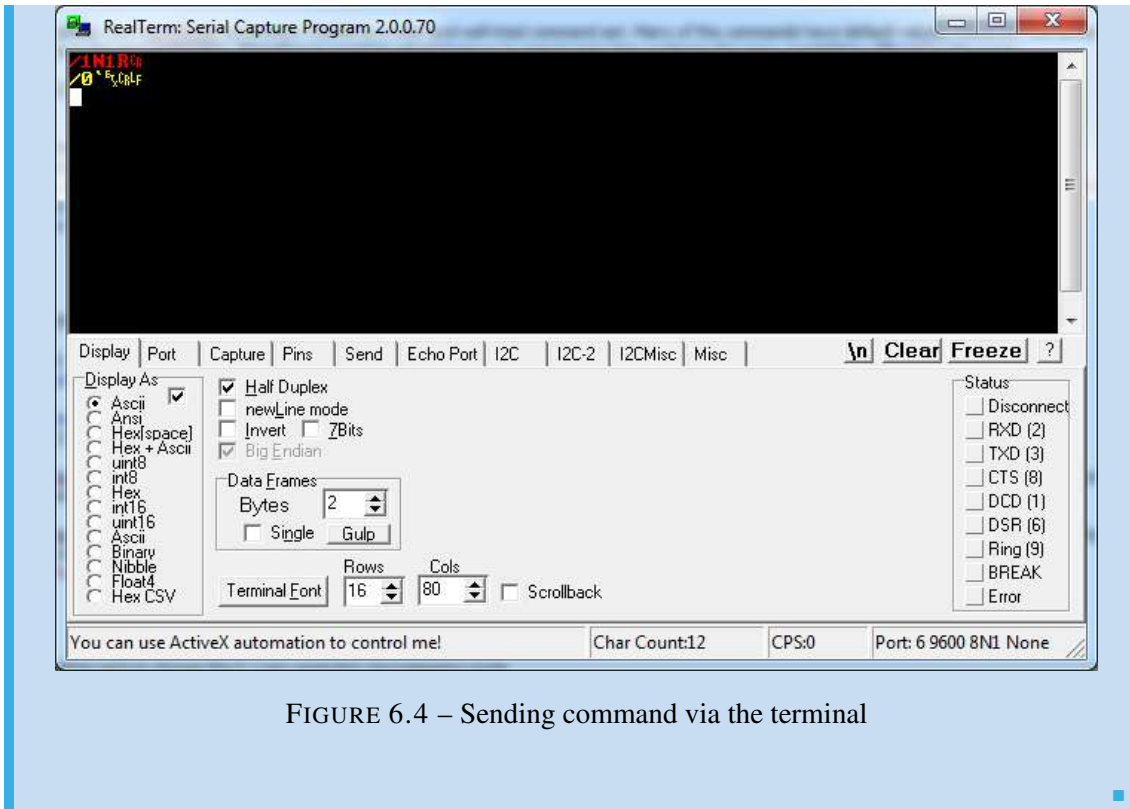
FIGURE 6.4 – Sending command via the terminal

*The speed setting is a delicate process since it is highly dependent on the microfluidics setup. Tubings of small inner diameter will act as restrictions for the flow and pushing liquids at high speed through such tubings will lead to high pressure in your microfluidic circuit.*

*High pressure will increase the leakage in the circuit and could even block the pump as too much force would be required to dispense the syringe content. Last but not least, high pressure can blow up the weakest part of a microfluidic setup. Therefore we recommend choosing speeds according to a pressure estimation.*

Speed Code	Value (pulses/sec)	Seconds/stroke (N=0, N=1)
10	1600	1.885
11	1400	2.15
12	1200	2.5
13	1000	3
14	800	3.75
15	600	5
16	400	7.5
17	200	15
18	190	15.79
19	180	16.665
20	170	17.645
21	160	18.75
22	150	20
23	140	21.43
24	130	23.075
25	120	25
26	110	27.275
27	100	30
28	90	33.335
29	80	37.5
30	70	42.855
31	60	50
32	50	60
33	40	75
34	30	100
35	20	150
36	18	166.665
37	16	187.5
38	14	224.235
39	12	250
40	10	300

TABLE 6.2 – Speed code for plunger movement

To better understand the relations between the motor speed, flow rate, plunger displacement and volume dispense, please refer to the appendix called “Resolution” in the user manual.

**Report Commands**

These commands do not need a trailing [R] character.

Command	Description
Q	Current status
? or ?0	Report absolute plunger position
?2	Report maximal speed (unit changes with speed mode)
?4	Report actual position of plunger
?5	Report speed mode (u = 0, U = 1, V = 2)
?6	Report valve position
?17	Report number of valve movements
?18 or %	Report number of valve movements (since last report)
?20 or #	Report firmware checksum
?23 or &	Report firmware version
?25	Report acceleration slope setting (plunger movement)
?26	Report pump address
?27	Report deceleration slope setting (plunger movement)
?28	Report current mode (fine positioning or microstep)
?29	Same as Q (query, status and error bytes)
?76	Report pump configuration
?300	Report plunger current (x10 mA)
?500	Report answer mode
?801	Number of valve positions
\$	Internal reset
*	Report supply voltage (x0.1 V)
?9000	Unique ID
?9010	Answers 1 if pump is initialized, 0 if not
?9100	Detailed status of the syringe
?9200	Detailed status of the valve

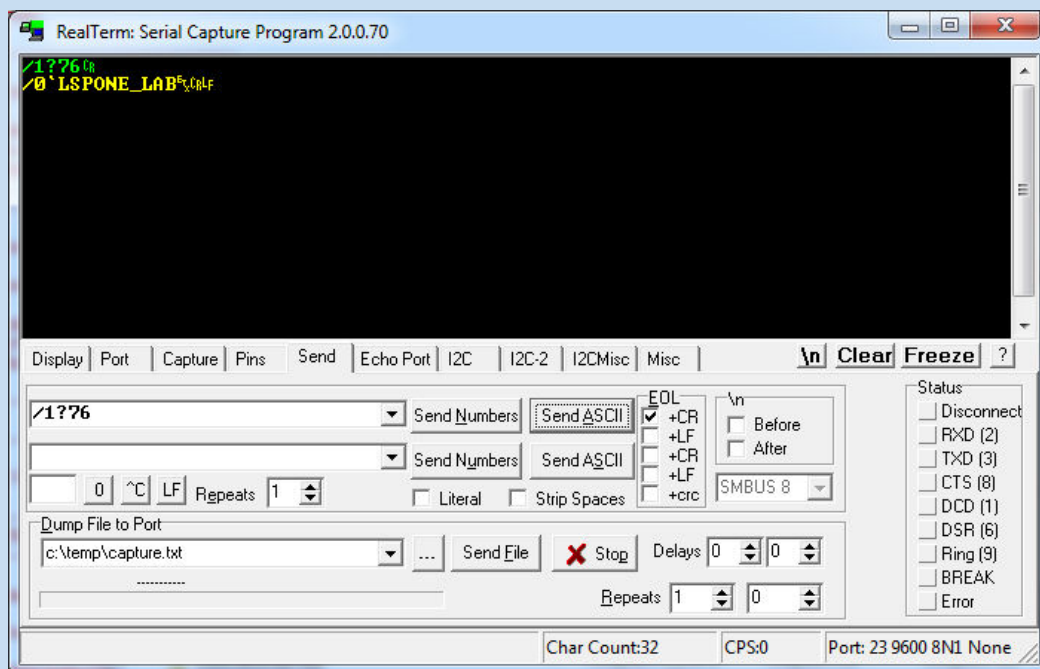
FIGURE 6.5 – Asking for the pump configuration

### Example 6.6

## 6.2 Debugging software problems

### 6.2.1 Visual feedback to the user

A very simple check when writing a sequence for the LSPone syringe pump can be obtained by looking at the Power-On button.

- When the blue LED is on, the pump is ready to receive commands.
- When the LED is blinking at 2 Hz, the pump is busy and the command required is being completed.
- When the LED is blinking at 5 Hz, an error has been encountered, the pump requires an initialization
- When the LED is blinking with an irregular pattern, an error has been encountered, the pump does not require an initialization

The fast blinking rate is the first step to debug any problem the user is facing when using the pump.

*Before looking in more detail at what caused an unexpected behavior of the pump, please make sure that:*

- *The pump is correctly powered. If the Power-On button emits no light at all, please check the power supply and the electrical connections.*
- *Check the hardware situation. The valve, the syringe and the plunger must be correctly tightened in their respective position.*
- *Make sure that you have sent an initialization command before running any other instructions.*
- *Do not forget to address the command to the pump by writing “/I<command>” (if the pump has the address “1”, in case you are not sure of the pump address, use “\_” which is the broadcast address) otherwise the pump will not be affected by the message.*
- *Do not forget to add R at the end of your command and to communication using the carriage return <CR> or \r*

After checking these first level errors, it is possible to explicit the error message using the report commands.

### 6.2.2 Error Codes and Pump Status

The pump reports immediately after it receives a command. This principle allows to check whether the command was accepted or not by the pump or whether it encountered any problem during the task.

**Note — Coding best practice.** A good coding practice when using this communication protocol is to check the pump status and catch possible errors with the [Q] command after each instruction.

#### Error Byte description

The error code has a length of 1 byte, i.e. 8 bits. The bit 5 is the status bit and serves for indicating whether the pump is busy or not. The error code is included in bits 0-3.

Bit	7	6	5	4	3	2	1	0
Value	0	1	Status Bit	0	Error Code			

TABLE 6.3 – Detail of the error byte as returned after a query “/IQ”

**Note — Interest of the (Q) query.** The answer block which is automatically sent by the pump features a status bit. However the status bit should not be used to know whether the pump is busy or not. A [Q] command is the only mean to know the current status of the pump.

### Status bit description

These statuses shown as a direct answer when you send a command to the pump.

Status Bit 5	Description
X=0	The pump will only accept report commands or terminate command [T]
X=1	The pump is ready to accept new instructions

TABLE 6.4 – Status bit description

**Note — Error byte after Terminate command.** After sending a terminate command, the status is busy but in this very case, the user can still send a new command

Error Byte 7 6 5 4 3 2 1 0	Symbol ASCII		Error	
	if Bit 5 = 0	if Bit 5 =1	Code	Description
0 1 X 0 0 0 0 0	@	‘	0	No Error
0 1 X 0 0 0 0 1	A	a	1	Initialization
0 1 X 0 0 0 1 0	B	b	2	Invalid command
0 1 X 0 0 0 1 1	C	c	3	Invalid operand
0 1 X 0 0 1 0 0	D	d	4	Missing trailing [R]
0 1 X 0 0 1 1 1	G	g	7	Device not initialized
0 1 X 0 1 0 0 0	H	h	8	Internal failure (valve)
0 1 X 0 1 0 0 1	I	i	9	Plunger overload
0 1 X 0 1 0 1 0	J	j	10	Valve overload
0 1 X 0 1 0 1 1	K	k	11	Plunger move not allowed
0 1 X 0 1 1 0 0	L	l	12	Internal failure (plunger)
0 1 X 0 1 1 1 0	N	n	14	A/D converter failure
0 1 X 0 1 1 1 1	O	o	15	Command overflow

TABLE 6.5 – Error codes

### Error Types

**Immediate errors** include Error 2 and 3. The automatic answer block sent after an instruction will feature the error character. No [Q] query is required to get a description of the error.

**Example 6.7** “/I014R<CR>” will trigger the following answer “/0c<ETX><CR><LF>” which means that an invalid operand has been used in the instruction. Here port “14” which does not exist on the valve. ■

To get detailed information about other errors, the user has to send the [Q] command and

analyze the subsequent answer. **Initialization errors** appear when the initialization could not be performed. Therefore the pump cannot be used until the error was cleared and a successful initialization completed.

**Command overflow errors** are caused by a Move command, a Set command (except [V]), or a Valve command being sent while the plunger is moving. The pump ignores the command and issues an error 15. The user should use the [Q] query to be informed when additional commands can be sent.

#### Error codes given by ?9100


Status code	Name	Description
FF = 255	Busy	Pump currently executing an instruction
00 = 0	Done	Pump available for next instruction
80 = 128	Unknown command	Check that the command is written properly.
90 = 144	Not homed	You forgot the homing! Otherwise, check that you have the right port configuration and try again.
91 = 145	Move out of range	You're probably trying to do a relative positioning and are too close to the limits.
92 = 146	Speed out of range	Check the speed that you're trying to go at.
E0 = 224	Blocked	Something prevented the plunger to move.
E1 = 225	Sensor error	Unable to read position sensor. This probably means that the cable is disconnected.

TABLE 6.6 – Detailed status of the pump (plunger) (see command ?9100)

## Error codes given by ?9200

Status code	Name	Description
FF = 255	Busy	Valve currently executing an instruction
00 = 0	Done	Valve available for next instruction
80 = 128	Unknown command	Check that the command is written properly.
90 = 144	Not homed	You forgot the homing! Otherwise, check that you have the right port configuration and try again.
E0 = 224	Blocked	Something prevented the valve to move.
E1 = 225	Sensor error	Unable to read position sensor. This probably means that the cable is disconnected.
E2 = 226	Missing main reference	Unable to find the valve's main reference magnet during homing. This can mean that a reference magnet of the valve is bad/missing or that the motor is blocked during homing.
E3 = 227	Missing reference	Unable to find a valve's reference magnet during homing. Please check that you have the correct valve number configuration with command "/1?801". If not, change it according to the valve you are working with. This can also mean that a reference magnet of the valve is bad/missing or that the motor is blocked during homing.
E4 = 228	Bad reference polarity	One of the magnets of the reference valve has a bad polarity. Please check that you have the correct valve number configuration with command "/1?801". If not, change it according to the valve you are working with. This can also mean that a reference magnet has been assembled in the wrong orientation in the valve.

TABLE 6.7 – Detailed status of the valve (see command ?9200)

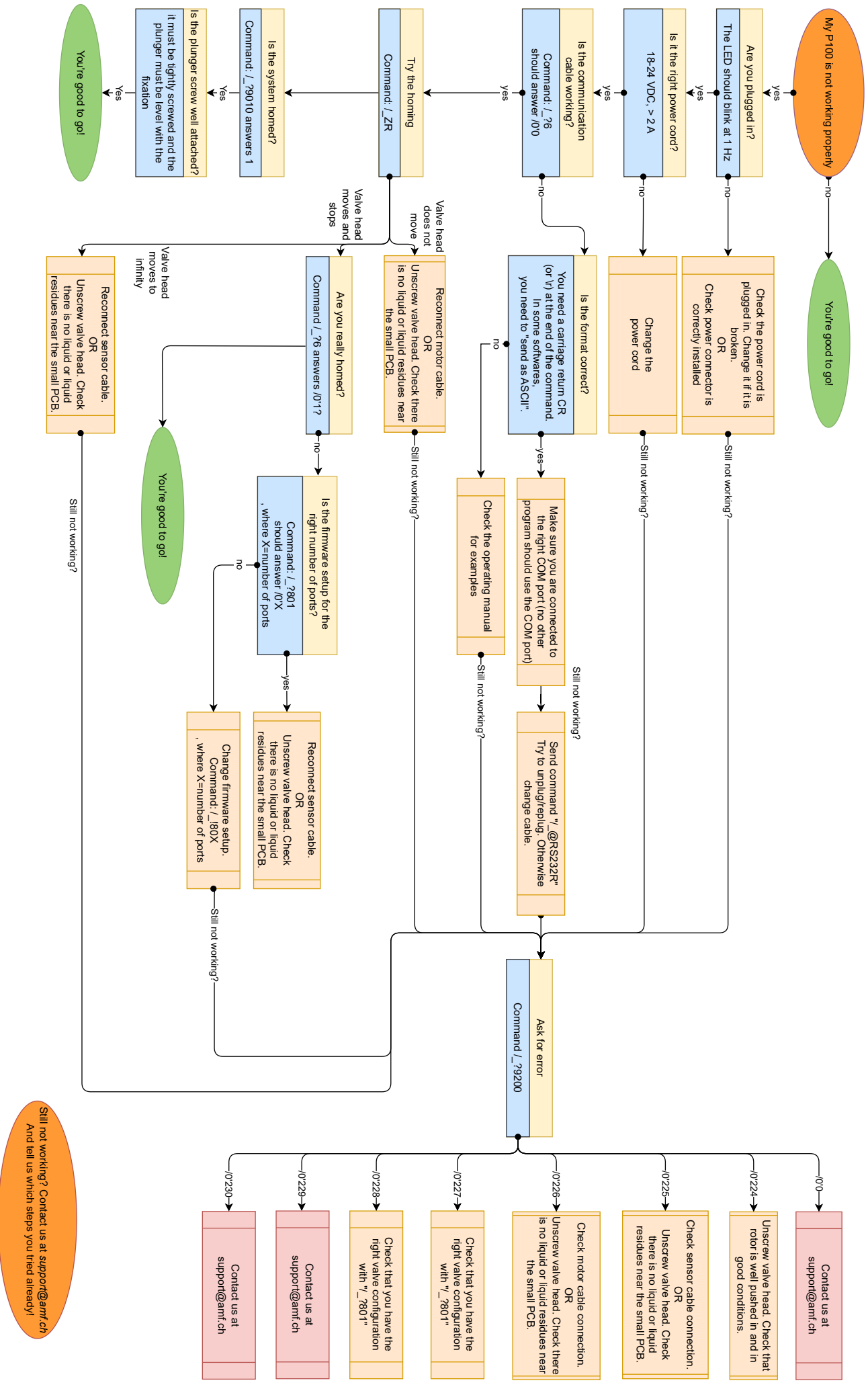## 7. GUI LSPoneQuick

### 7.1 Description

Offering an easy software tool to start experimenting with the LSPone syringe pump within a few seconds is a main commitment at *Advanced Microfluidics SA*. A cross-platform graphical user interface has thus been developed for a plug-and-play experience. Among other features, the program offers an automatic detection of the hardware as well as a macro recording tab to automate your most requiring sequences. Moreover, firmware updates can also be done within this environment. We can thus help our clients keep their device up-to-date without displacing their hardware.


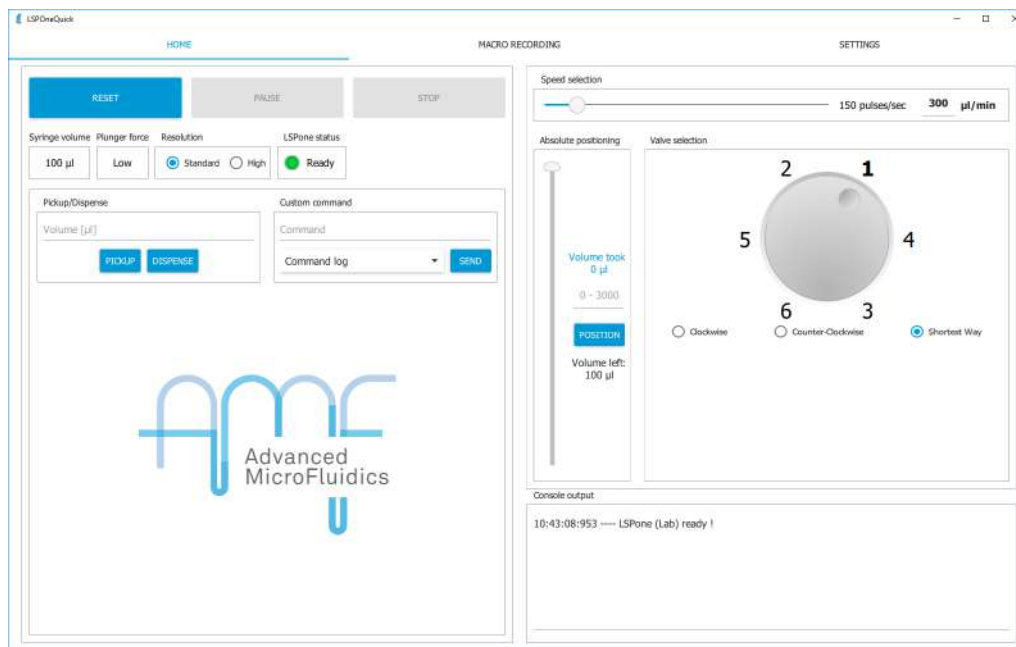
FIGURE 7.1 – Main panel of the Graphical User Interface


### 7.2 Requirements

The LSPoneQuick software is cross-platform and will thus work on most operating systems. The software has been successfully tested on Windows 7, 8 & 10 and Mac OS X.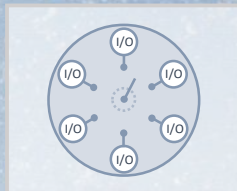

## 8. Labview VI
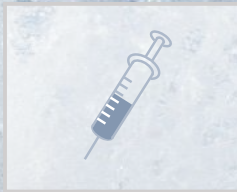
### 8.1 Description

For ease of use, a sub-VI in Labview was created. You can directly integrate it in your current Labview setup, providing the COM port and string command as input, and observe the response from the pump to check that everything is acting as wanted.

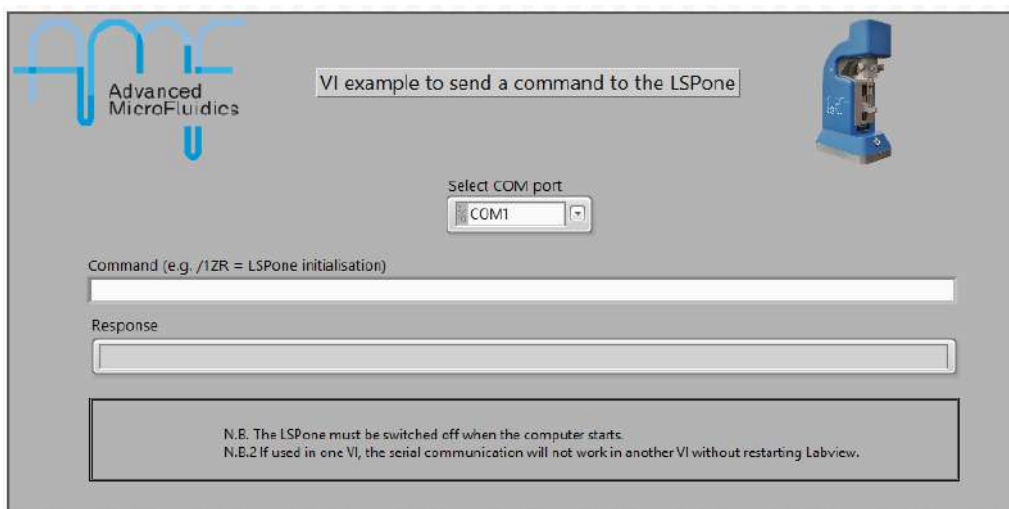
FIGURE 8.1 – Labview sub-VI, front panel - 1 pump

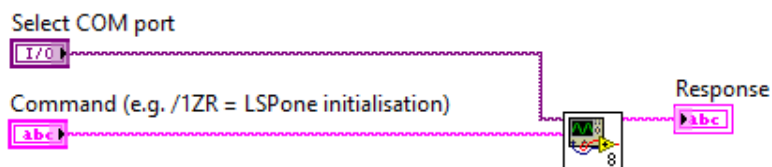
FIGURE 8.2 – Labview sub-VI, bloc diagram - 1 pump

It is then very simple to include several pumps in the same setup. In the following example, the same command is sent to 2 LSPones, but it is also possible to send 2 different commands.

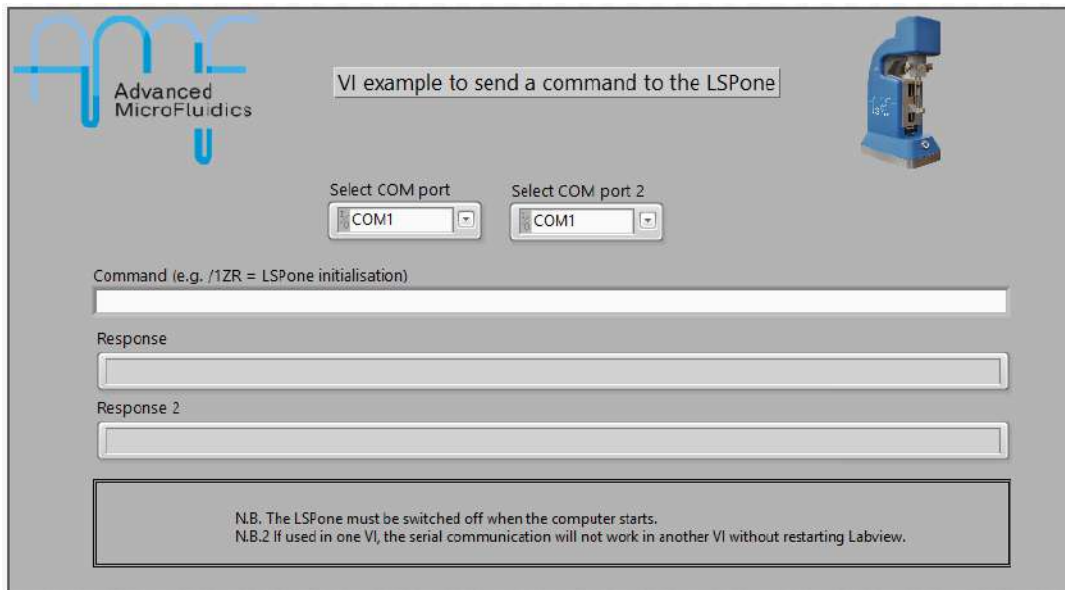
FIGURE 8.3 – Labview sub-VI, front panel - 2 pumps


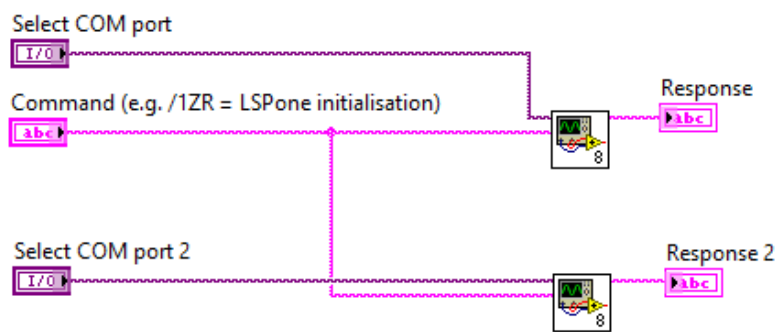
FIGURE 8.4 – Labview sub-VI, bloc diagram - 2 pumps


## 8.2 Requirements

Nothing special is required, other than Labview from National Instruments. Tested with Labview 2016.

*N.B. In case the VI does not seem to respond, please restart Labview.*

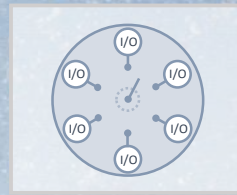## 9. Python scripting
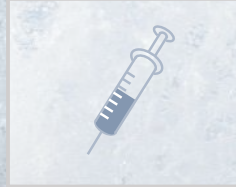
### 9.1 Description

Serial commands can be sent via python scripting using the *pySerial* library. A very simple example is shown below. Python scripting will be of specific interest when using different devices and trying to automate a sequence involving these devices.

```
#!/usr/bin/env python
'''
~~~~~
Script for LSPone - Example code
~~~~~
'''

import serial
import time

COM_PORT = "COM10" # WILL BE DIFFERENT ON YOUR SYSTEM
BAUDRATE = 9600
TIMEOUT = 1000 # [ms]

# # LSPone initialization
lsp = serial.Serial(COM_PORT, BAUDRATE, timeout=TIMEOUT)
print(lsp.name)
lsp.write(b"/1ZR\r")
```

FIGURE 9.1 – Python example script

### 9.2 Requirements

Python has to be installed on your computer. Additionally, the script requires the *pySerial* library (<https://pythonhosted.org/pyserial/>).

# A. Resolution

## A.1 Definitions

**Full stroke** Movement of the plunger from top to bottom (i.e. 30 mm)

**Pulse** Displacement of 0.01 mm of the plunger

**Step** Subdivision of a pulse

**Step resolution** Displacement of a step

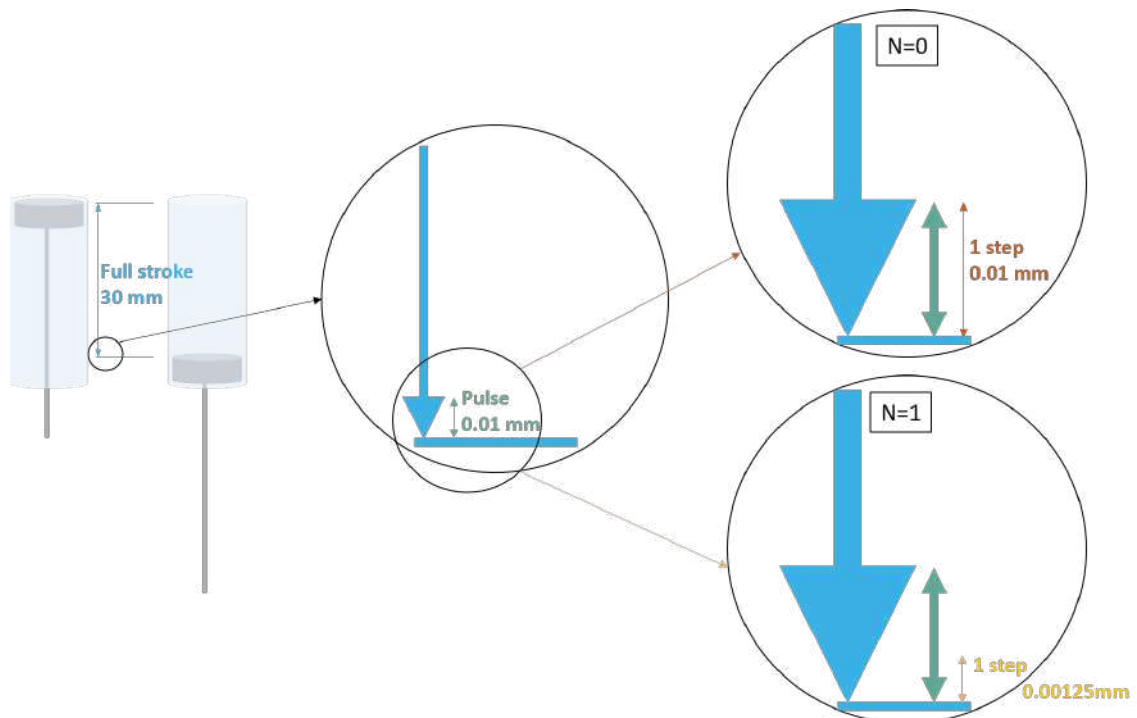
FIGURE A.1 – Definitions in image

## A.2 Flow rate in function of plunger speed

The speed of the plunger is always defined in pulses/s = displacement of the plunger per second. A relation can be found with the flow rate of the liquid.

Plunger speed [pulses/s]	Syringe size [ $\mu\text{L}$ ]						Flow speed [ $\mu\text{L}/\text{min}$ ]
	25	50	100	250	500	1'000	
0.00745 (speed "u1")	0.003725	0.00745	0.0149	0.03725	0.0745	0.149	
0.05 (speed "U1")	0.025	0.05	0.1	0.25	0.5	1	
5 (speed "V5")	2.5	5	10	25	50	100	
10	5	10	20	50	100	200	
50	25	50	100	250	500	1'000	
100	50	100	200	500	1'000	2'000	
500	250	500	1'000	2'500	5'000	10'000	
1'000	500	1'000	2'000	5'000	10'000	20'000	
1'500	750	1'500	3'000	7'500	15'000	30'000	

TABLE A.1 – Flow rate in function of plunger speed

**Example A.1** For a 100  $\mu\text{L}$  syringe

5 pulses/s = 10  $\mu\text{L}/\text{min}$  = 10  $\mu\text{L}/60\text{s}$  = 1/6  $\mu\text{L}/\text{s}$

**In 1 second, the plunger does 5 pulses (= 5\*0.01 mm), during which 0.17  $\mu\text{L}$  are delivered.**

■

## A.3 Volume in function of plunger displacement

The position at which the plunger can stop can be changed, i.e. a pulse can be subdivided into steps of various sizes. This allows us to attain different precisions in the quantity of volume delivered.

Resolution type	Resolution command N	Resolution [steps/stroke]	Resolution [steps/pulse]	Resolution [mm/step]
Standard	N=0	3'000	1	0.01
High	N=1	24'000	8	0.00125

TABLE A.2 – Resolution in different units

Resolution [steps]	Syringe size [ $\mu\text{L}$ ]						Volume delivered for 1 step [nL]
	25	50	100	250	500	1'000	
3'000	8.33	16.67	33	83.3	166.7	333	
24'000	1.04	2.08	4.2	10.4	20.8	41.7	

TABLE A.3 – Minimal volume delivered in function of plunger course resolution

**Example A.2** For a 100  $\mu\text{L}$  syringe:

100  $\mu\text{L}$  = 3'000 steps => 1 step = 0.033  $\mu\text{L}$  = 33 nL

100  $\mu\text{L}$  = 24'000 steps => 1 step = 0.0042  $\mu\text{L}$  = 4.2 nL

**In standard resolution, 1 step corresponds to a volume of 33 nL.**

**In high resolution, 1 step corresponds to a volume of 4.2 nL.**

■