

SPM HD/P110-O Programmable Syringe Pump

Low Volume - DATASHEET




Advanced Microfluidics

Environmental conditions

Operating temperature	15-45°C (59-113°F)
Humidity	20-80%, non-condensing
Max pressure	7 bars (102 psi)

Fluidic characteristics

Tube port fittings	Standard ¼-28 UNF, flat-bottom
Wetted materials	PTFE or UHMW-PE, PCTFE and borosilicate glass
Channel diameter	0.5 mm (0.020 in) / 0.4-1 mm (0.015-0.039 in) available upon request

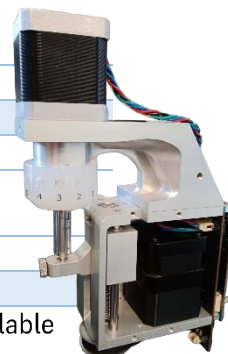

Mechanical characteristics

COMPATIBLE VALVE MODELS

Series	ULTRA-LOW CARRYOVER VOLUME	LOW CARRYOVER VOLUME
Liquid path		
Number of channels inside rotary part	2	1
Number of ports	6	6, 8, 10, 12
Channel diameter	0.5 mm	0.5, 1 mm
Internal volume	5.2 µL	3.5 – 14.9 µL
Carryover volume	1.5 µL	2.6 – 7.8 µL
Dead volume	None	

Glossary reminder

Internal volume: Volume inside the system, from entrance to exit (does not include the volume in the syringe)

Dead volume: Volume that is “stuck” in the system (dead end), which is not cleanly swept and relies on diffusion to clear out.

Carryover volume: Volume of liquid that will be mixed with the next liquid or sample. It is not stuck, but will be swept next time a liquid passes.

PLUNGER DETAILS

Travel length	30 mm with 1'296'000 micro-steps
Resolution	Selectable 3'000 steps or 24'000 steps
Drive	Screw drive with linear encoder for step loss detection
Plunger seal material	PTFE (or UHMW-PE upon request)
Accuracy	≤ 1.0 % at full stroke

Advanced Microfluidics SA
Chemin de la Dent d'Oche 1a
CH-1024 Ecublens, Switzerland
Phone: +41 (0)21 552 14 30
Email: info@amf.ch
Web: www.amf.ch

SPM HD/P110-O Programmable Syringe Pump

Low Volume - DATASHEET




Advanced Microfluidics

COMPATIBLE SYRINGE MODELS

Ref. #	Plunger material	Syringe volume [μ L]	Min. flow rate [μ L/min]	Max. flow rate [μ L/min]	Min. dosing volume [μ L]
S-50-P	PTFE	50	0.00717	400	0.1
S-100-P	PTFE	100	0.0144	800	0.2
S-100-U	UHMW-PE	100	0.0144	800	0.2
S-250-P	PTFE	250	0.0359	2'000	0.5
S-500-P	PTFE	500	0.0717	4'000	1
S-500-U	UHMW-PE	500	0.0717	4'000	1
S-1000-P	PTFE	1'000	0.143	8'000	2

COMPLETE MODULE (Ref. P110-O)

Dimensions	199.7 x 136.7 x 49.6 mm
Weight	1.5 kg
Default configuration	500- μ L syringe with PTFE plunger (S-500-P) 6-position valve with low carryover volume (V-D-1-6-050)
External fixation system	4x M3 screws, position shown in orange

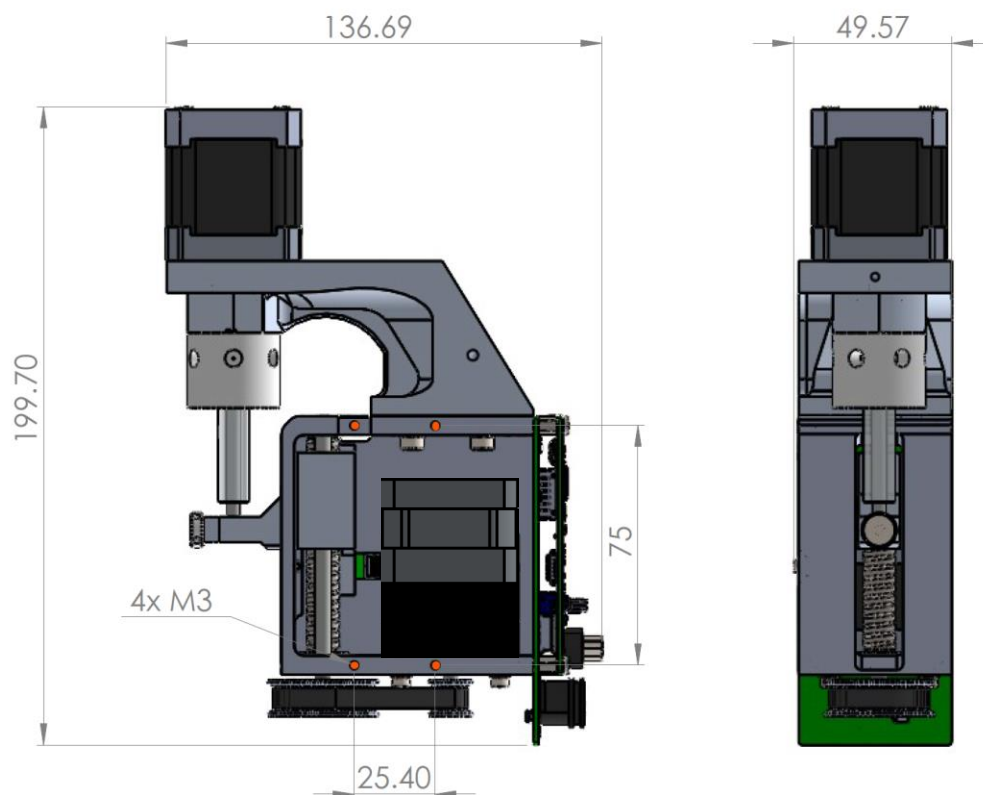

Dimensions are in mm

Electrical characteristics

Power	18-24 VDC, 2.2 A peak
Required cables for operation	USB mini to USB, Power cord (power jack 2mm)

Advanced Microfluidics SA
Chemin de la Dent d'Oche 1a
CH-1024 Ecublens, Switzerland
Phone: +41 (0)21 552 14 30
Email: info@amf.ch
Web: www.amf.ch

SPM HD/P110-O Programmable Syringe Pump

Low Volume - DATASHEET



Communication interface

Electrical Interface	USB mini, 9-pin D-Sub (other upon request)
Interface	USB mini, RS-232 or RS-485
Communication type	Serial communication

The valve controller is seen as a virtual serial port. There is an on-board FTDI USB to serial chip, so the drivers are automatically installed on Windows 7 and above.

SERIAL COMMUNICATION PARAMETERS

Baud rate	9600
Data Bit	8
Parity	None
Stop Bit	1
Handshake	None
End line	<CR>

View the SPM operating manual for the extensive command set.

Revisions

Version	Author	Release Date	Modifications	Revised By	Approved By
01.01	RRY	15/05/2025	First release	CAL	RRY
01.02	RRY	19/05/2025	Title update	NFA	RRY
01.03	RRY	26/05/2025	Update volume	MGE	RRY
01.04	MGE	15/07/2025	Update flow rates	CAL	MGE

The information and data contained in this documentation is subject to change without notice. The reproduction, distribution, and utilization of this document as well as the communication of its contents to others without explicit authorization is prohibited.

All rights reserved in the event of the grant of a patent, utility model or design.

The general terms and conditions of Advanced Microfluidics SA apply. Alternative agreements must be in written form.

Advanced Microfluidics SA shall not be liable for errors contained in this document or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.